



Technology for Productivity

Spectra™ Products

EasyView & e-Vision

As part of our continuous product improvement policy, we are always pleased to receive your comments and suggestions about how we should develop our product range. We believe that the manual is an important part of the product and would welcome your feedback, particularly relating to any omissions or inaccuracies you may discover.

You can send your comments to:-

MKS Instruments, Spectra Products
Cowley Way
Crewe
Cheshire
CW1 6AG
U.K.

+44 (0) 1270 250150 Tel.

+44 (0) 1270 251939 Fax.

Email – manual_info@mksinst.com
www.mksinst.com

e-Vision®, e-Vision+® and EasyView® are registered trademarks of MKS Instruments
Windows and Windows 95 are trademarks of the Microsoft Corporation.

All other brand or product names are trademarks or registered trademarks of their
respective companies.

<i>i. Getting Help</i>	5
<i>Warning</i>	6
Connecting Control Units.....	7
2. Installation Requirements	8
3. Installation	9
Overview.....	9
3.1 Installing EasyView	10
EasyView Set-up dialog.....	10
On-line dialog.....	12
Destination Folder dialog.....	12
Support Information dialog.....	13
3.2 Creating User Accounts	14
Overview.....	14
3.3 Creating the "RGA User" Account.....	15
4. The Set-up Wizard	20
Overview.....	20
4.1 Standard setup.....	21
EasyView Setup Wizard dialog.....	22
Hardware Setup dialog.....	23
Pressure Units dialog.....	24
Disk Storage Location dialog.....	24
Log File Details dialog.....	25
5. Adding your Instruments	27
5.1 Recommended installation method – If 192.168.0.250 is not in use.....	28
5.2 Recommended installation method – If 192.168.0.250 is already in use.....	29
5.3 Device Manager.....	30
Information pane.....	30
Controls.....	30
IP Configuration dialog.....	32
Device Manager Controls.....	33
Access Levels.....	35
6. Using EasyView	37
Overview.....	37
6.1 Starting EasyView.....	38
6.2 Toolbars.....	39
6.3 Analog Mode.....	41
6.4 Barchart Mode.....	43
6.5 Leak Check Mode.....	45
6.6 Peak Jump Mode.....	47
6.7 Disk Store.....	50
6.8 The Main EasyView screen.....	52
1 Title Bar.....	52
2 Menu Bar.....	52
3 EasyView Control Icons.....	52
4 Views Windows.....	52
5 EasyView Control Window.....	53

6 Toolbars	53
7 Default Toolbar	53
8 Scan Navigation Toolbar	54
9 Chart Display Toolbar	55
10 Trend Display Toolbar	56
Customizing Toolbars	56
7. Changing/Adding Views in EasyView	61
7.1 Swept Measurement Properties Settings	63
X-Axis.....	63
Initial Y-axis Properties	64
7.2 Trend View Properties Settings	66
Y-Axis Auto-range Properties.....	74
8. Other Menus	83
View Pull-Down Menu	83
Toolbars	83
Status Bar.....	83
Properties	84
Window pull-down menu	84
New Trend Window	84
Arrange All Windows	85
Bring Recipe to Front.....	85
Cascade Recipe	85
Tile Recipe Horizontally.....	86
Tile Recipe Vertically	86
Selecting Individual View	86
9. EasyView Diagnostics	87
10. EasyView Tuning	89
11. EasyView Calibration	91
12. Preferences.....	98
Units of Pressure.....	98
Auto View to front.....	99
Graph Settings	99
Common Gases List.....	99
13. Printing.....	100
Printing a View	100
14. EasyView Tools.....	103
Backup and Recover	103
Computer Lock	106
Exiting Computer Lock.....	108
15. Java Web Applet.....	110
15.1. Accessing the instrument.....	111
15.2 Page Views and Controls	114
The Tool Bar	114
Page Tabs.....	116


i. Getting Help

We are always pleased to provide assistance where we can. If you are experiencing difficulties or need help, feel free to call your local MKS Instruments facility and ask for the Customer Support Group.

Please have the following information ready so that our technical staff may help you quickly and effectively:

The serial numbers of the analyser and e-Vision unit.

The EasyView software version number, this is written on the software CD ROM and can also be found when using EasyView by clicking **Help/About**.

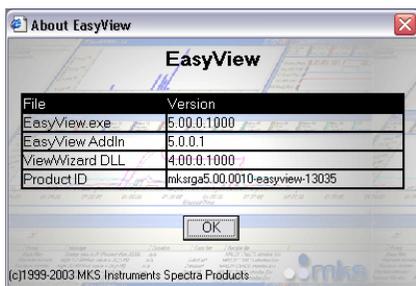

The operating system used.

On-line help is available during the install process and can be accessed by clicking the <Help> button on any of the dialog boxes.

Support Contact Telephone Numbers:

United Kingdom
+44 (0) 1270 250150

Warning

The control unit must not be powered up when not connected to the analyser

1. Introduction

EasyView is designed to operate with the MKS Spectra e-Vision range of instruments. It will not operate with any other control unit in the MKS Spectra range.

EasyView will run under Windows 98SE/ME/NT/2000 and Windows XP operating systems. No other versions of Windows are supported.

The EasyView package is designed to monitor partial pressures in a production environment where it is essential to observe and record the current status of the process, while simultaneously reviewing historical events in both the current and archived runs.

The data acquisition is fast and covers a wide dynamic range of pressures so that some or all of the partial pressures can be measured and stored, while only data of interest is displayed.

The package has been designed to be easy to use and understand so it can be used by skilled and unskilled operators alike

Connecting Control Units

The e-Vision range uses an RJ-45 Ethernet connection only. The cable used must meet Cat 5e specification and be of the STP (Shielded Twisted Pair) type.

For connection of an e-Vision unit directly to the PC, the cable must be a cross-over type cable.

If the connection is made through a hub or network, the cable is a standard patch cable.

The control unit must be one suitably configured to run with EasyView. If the control unit was supplied with another software package or a previous version of EasyView, please contact your local MKS Spectra facility for upgrade instructions.

Note: This manual is concerned ONLY with the EasyView software package used in conjunction with the e-Vision range of instruments.

2. Installation Requirements

EasyView operates under Windows 98SE/ME/NT (SP6 or higher)/2000 and Windows XP environments, running on an IBM compatible PC.

It is recommended that ALL available updates for your particular operating system have been applied before the installation of EasyView.

The EasyView Installer will search your system for files required for the correct operation of the program, these may include Service Packs, Windows Installer Updates or newer versions of Internet Explorer.

This process is automated and simply requires the user's permission to install the required components.

Note: You will need to have administrative privileges to install these updates.

Once the required updates have been applied, or if your machine already meets the software requirements, the first page of the EasyView Installation Wizard appears.



See Section 3 for details on installing EasyView.

3. Installation


Overview

Before you start to install EasyView, make sure you have fitted and configured any network cards you may require.

If you are using Windows NT/2000 or XP, you must log on as an Administrator to install EasyView. You will only require "Power User" privileges to run EasyView.

This manual describes each stage of installation in turn and each section should be completed fully, before moving onto the next.

Enter a "Full Name" and "Organization" if required, otherwise leave at the default.

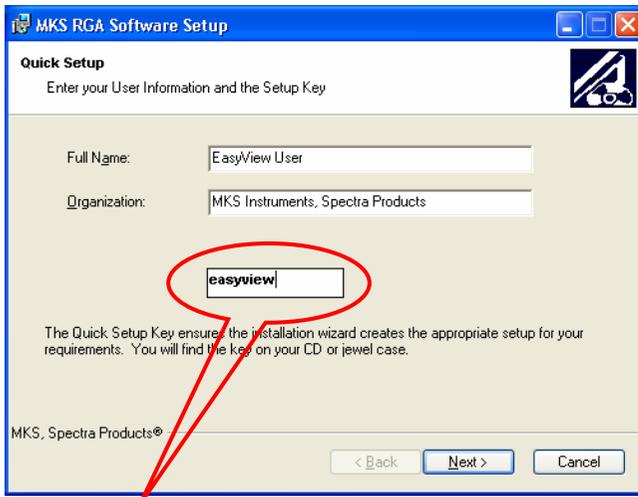


The "Quick Setup Key" is supplied with your CD and determines the type of installation.

For an EasyView installation, the CD key is "easyview".

If an invalid key is entered, a confirmation dialog appears giving you the option to re-enter the key.

Once a valid key has been entered, click the <Next> button to continue.


On-line dialog

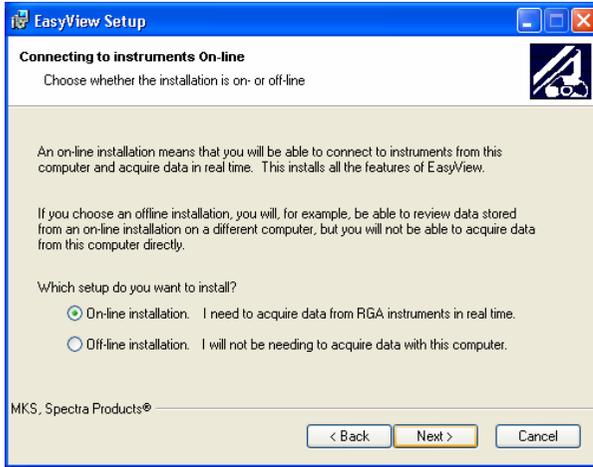

An “On-line” installation installs the full suite of EasyView components and should be chosen if the PC is to run the RGA instrument

An “Off-line” installation installs only the components required to view stored data and create or modify recipes.

Destination Folder dialog

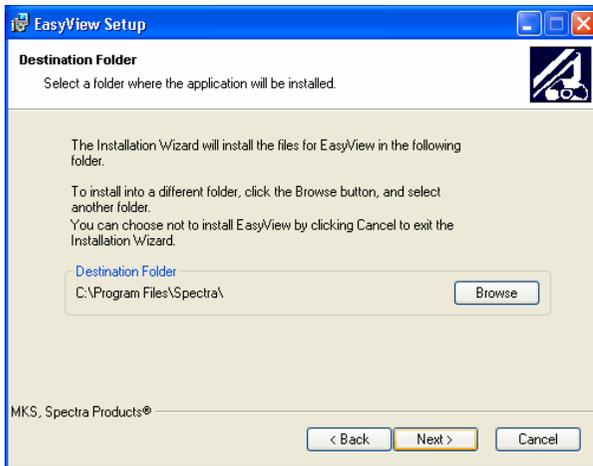


Here you can select the installation directory used for the EasyView software. The above shows the default directory of **C:\Program Files\Spectra**.

If you wish to install to another directory, click on the <Browse> button. A standard "Browse" dialog is displayed where you can select a directory of your choosing.

The installation process will begin to install the EasyView software onto your PC. This process may take a few minutes, depending on the type of installation chosen and the specifications of the PC.

Support Information dialog

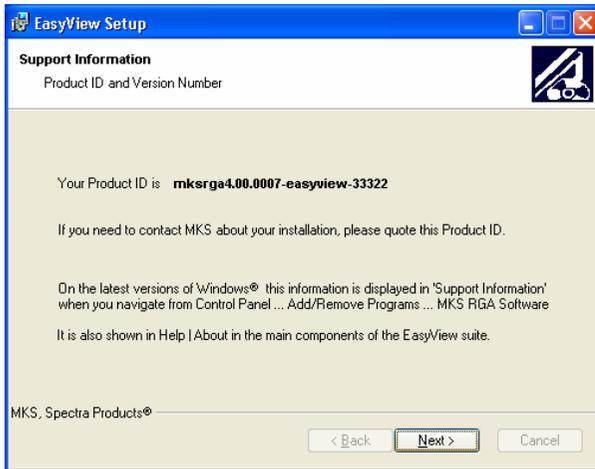

The "Support Information" dialog provides details of your software version numbers and instructions on how to re-obtain these version numbers after installation.

Note: the version number will be required if support is requested.



Once the program is installed you should work through Section 3.2 Creating User Accounts.

3.2 Creating User Accounts

Overview


The following section describes how to create a new user and assign user rights using the WindowsXP environment as a guide.

This is not common for all Windows operating systems, advice from your IT representative or MKS Spectra should be sought if you are unsure of how to continue, or if the PC is part of your company network.


You will need to create a "RGA User" account with "Power User" rights.

3.3 Creating the “RGA User” Account


Click on the <Start> button, right-click on “My Computer” and select “Manage” from the menu.

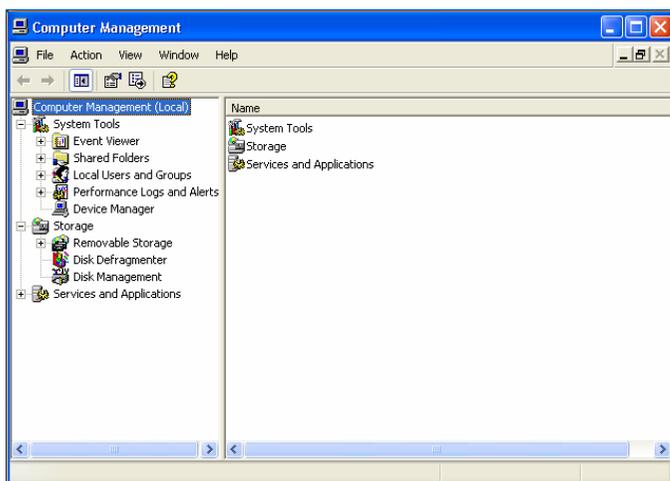

In the “Computer Management” dialog, click on the **Local Users and Groups** entry to expand the list and show the **Users** and **Groups** folders.

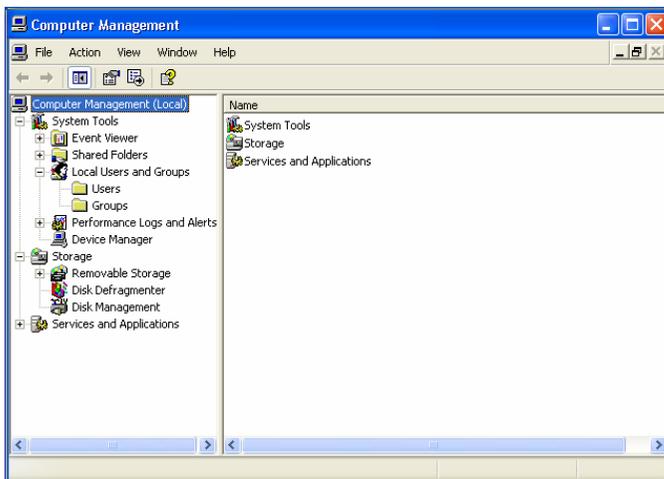


Right-click the **Users** folder and select **New User** from the list.

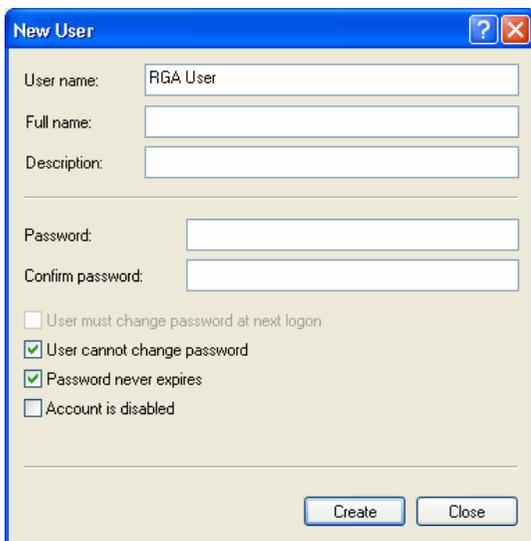


It is recommended that you use the name "RGA User" to easily identify the account. You can enter a full name, description and password if you wish.

Check the **User cannot change password** and **Password never expires** boxes as shown.


Click on the <Create> button and then click the <Close> button to return to the "Computer Management" dialog.

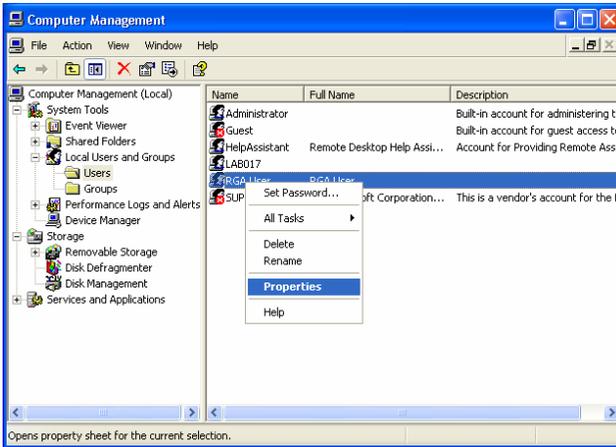


Right-click the "RGU User" account you have just created and select **Properties** from the list.

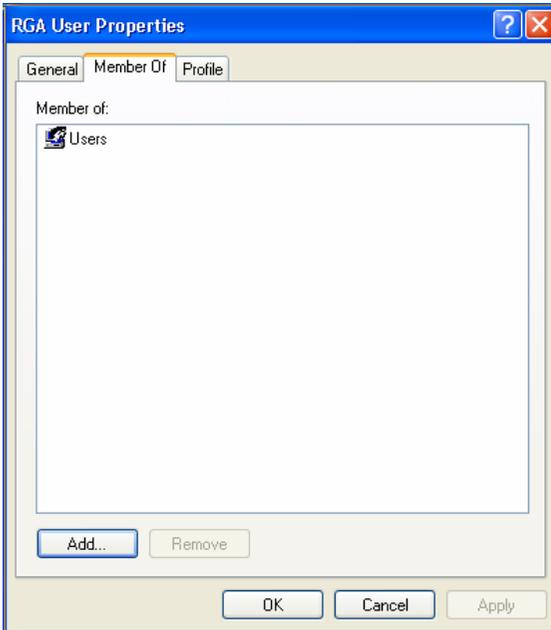


Click on the **Member of** tab to display the default **Users** group. Click on the <Add> button to add another group type.

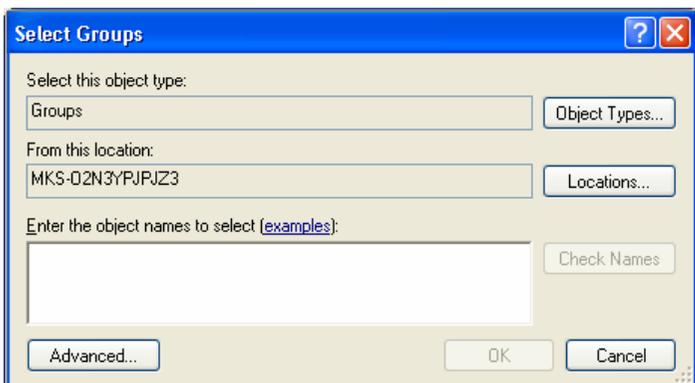


In the **Enter the object names to select** field, type in "Power Users" as shown and click <OK>.

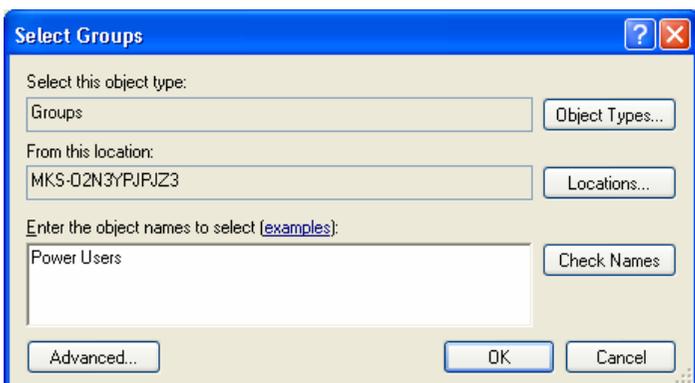

The **Select Groups** dialog will reflect the changes made.

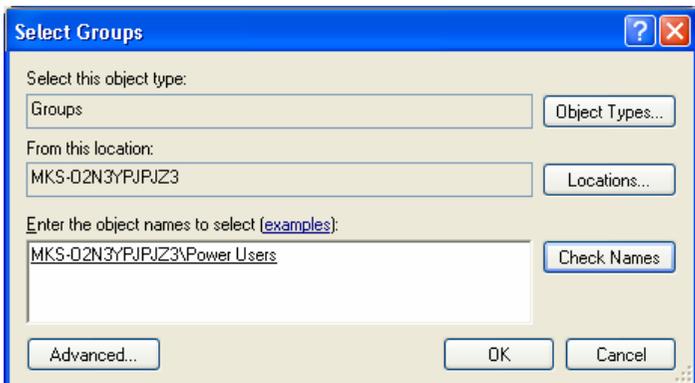


Click <OK> to return to the “Computer Management” dialog.

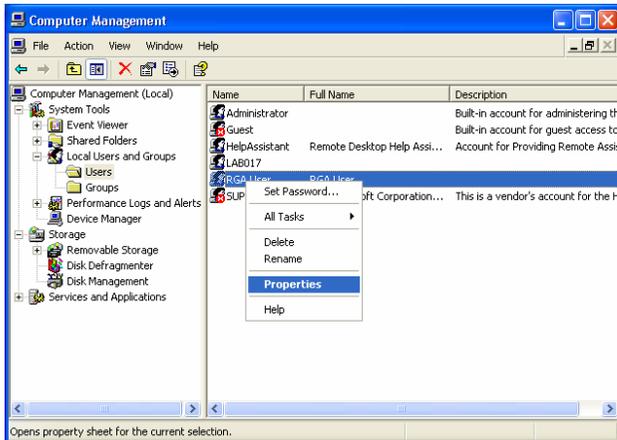

Right-click on the **RGU User** account and select the **Properties** field.

In the “RGU User Properties” dialog, confirm that **Power User** is displayed under the **Member of** tab.

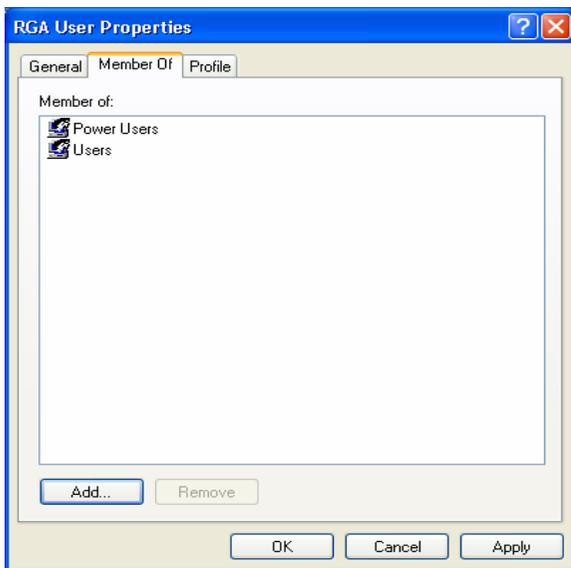

Click on the <OK> and exit the “Computer Management” dialog.

At this point, you should now log-off as the Administrator and log-on as the RGA User to continue through Section 4. - Set-up Wizard.

4. The Set-up Wizard

Overview


The Setup Wizard application configures the connection to the instrument, the directory used to store data, default pressure units and how the EasyView software will start.

Once EasyView has been installed and the RGA User account created, you must run the "Setup Wizard" before you can use the software.

This can be found in **Start/All Programs/MKS RGA Applications/Tools/Setup Wizard**, or by attempting to start the EasyView program prior to running Setup Wizard.

4.1 Standard setup

Click on **Start/All Programs/MKS RGA Applications/Tools** and select the **Setup Wizard** entry.

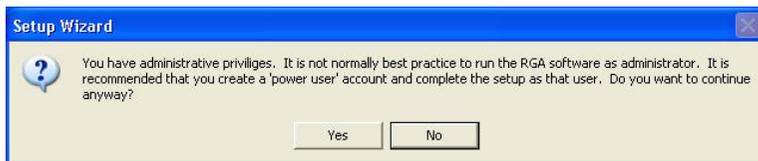

If you are currently logged on as an administrator, the above warning dialog will appear. It is best to log-off and switch accounts to the RGA User you created earlier.


If an earlier version of our RGA software is detected by the Setup Wizard, the following dialog is displayed.



This manual is concerned only with e-Vision instruments. Only the option shown above should be selected.

Do not attempt to carry out any upgrade to any existing non e-Vision based instruments you may have.


EasyView Setup Wizard dialog

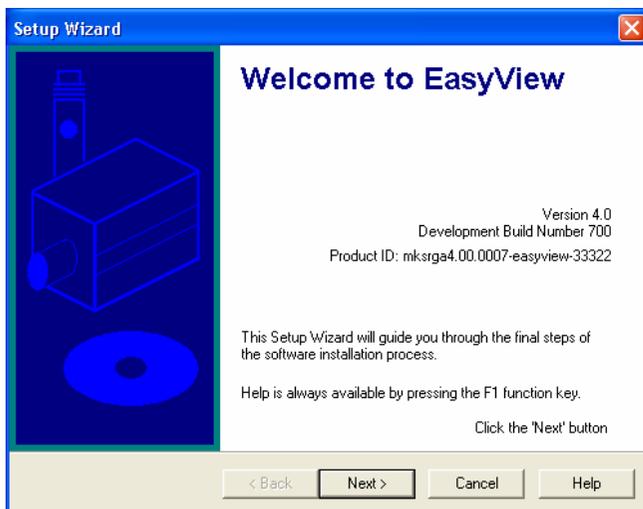

At any stage during software setup you can click on <F1> to bring up the help dialog.


These help pages give concise information about the current page of the setup wizard.

Hardware Setup dialog

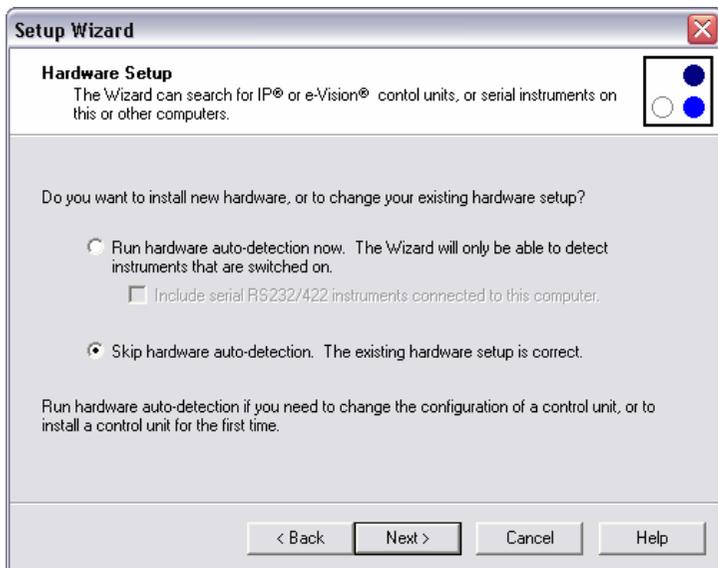

For now, click the option to skip hardware auto-detection as shown and continue to setup the EasyView software.


Pressure Units dialog.

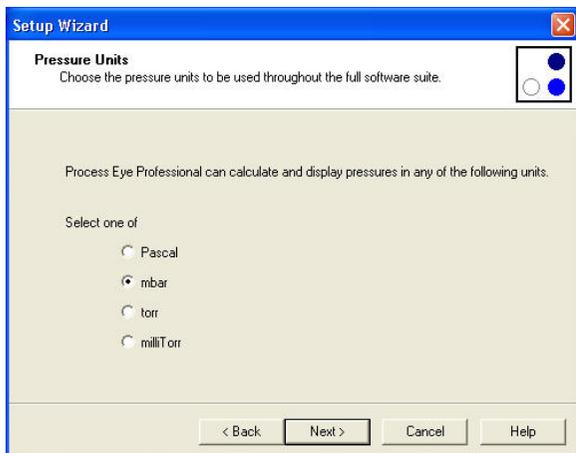

Here you may choose the unit of pressure EasyView will use by default.

This may also be changed later while using EasyView, by selecting **Recipe / Preferences** on the menu bar.

Disk Storage Location dialog.

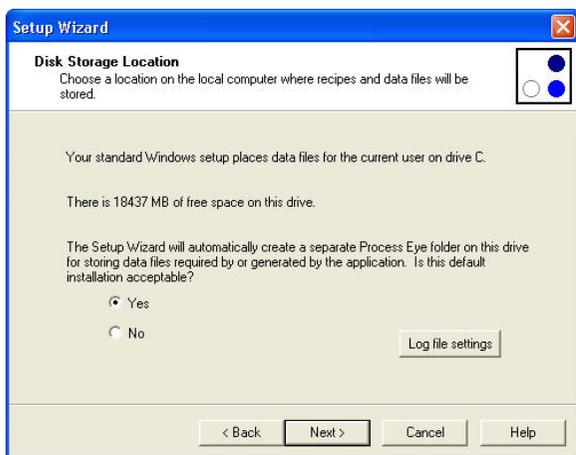

The default location for storing recipes, views and data is **Current User/My Documents**.

If multiple users are to use the same PC, logging on with different usernames, it may be simpler to create a common location available to all users.

Choose <No> to specify a folder of your choosing.


Log File Details dialog.

The "EasyView Error Log File" is where information on any script errors encountered is stored. This information can be viewed to help with recipe diagnosis.

The default value is adequate in most cases, but if you need to increase the file size, or store the log as an XML file, these options can be set here.



Otherwise, accept the default recommendation and click <Next>.

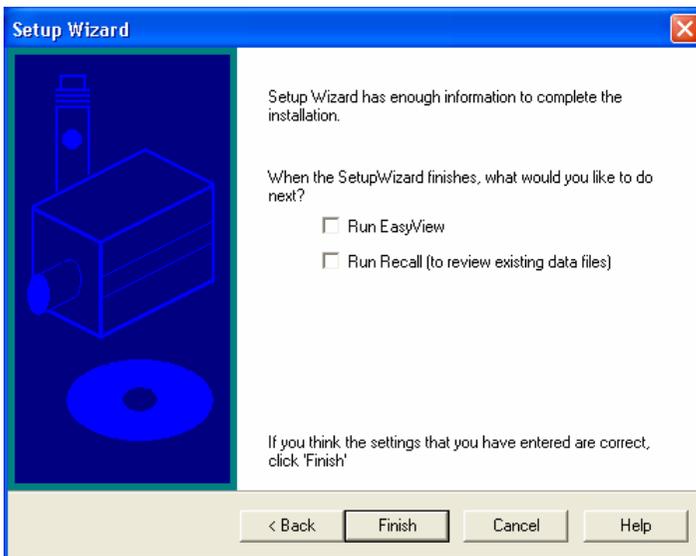

To complete the standard installation, leave the boxes un-checked and click the <Finish> button.



An "installation completed" dialog will appear, click <OK> to exit the installation program.

5. Adding your Instruments

Before connecting any instruments to your network, you must seek advice from your IT staff on the following points:

Is the IP address 192.168.0.250 already in use?

All new instruments have a default, fixed IP address of 192.168.0.250 which can be changed to suit your network if required. Only one instance of an IP address can exist at any one time, this address needs to be available for the initial configuration of the instrument.

What IP addresses are available for my instruments?

You will require a unique IP address for each instrument.

What is the Subnet range of my network?

This will be required during the configuration of the instrument.

Concise help on IP addresses, networking, network configuration and using Device Manager can be accessed by either choosing the <Help> button on the Device Manager dialog, or from the MKS RGA Applications program group.

We recommend you read through these help pages before continuing any further.

5.1 Recommended installation method – If 192.168.0.250 is not in use

After obtaining the information you required above, you should connect your instruments to your network using the supplied Cat5e shielded UTP cable. If this cable is not a suitable length, an alternative of the same specification should be sought.

Assign a new IP address to each instrument in turn and then install the instruments on your network.

Proceed to Section 5.3 - Device Manager.

5.2 Recommended installation method – If 192.168.0.250 is already in use

If the above IP address is not available on your network, then the instrument must be configured using a PC disconnected from the network.

Seek advice from your IT staff on obtaining a PC with a fixed IP address within the range of addresses used by your network, but not one that you intend to use for an instrument.

Connect the e-Vision directly to the PC using a “Cross-over” cable, or if you do not have one of these cables available, use a standard patch cable and a hub.


Assign a new IP address to each instrument in turn and then install the instruments on your network.

Proceed to Section 5.3 - Device Manager.

5.3 Device Manager

Before you can begin to use an instrument with EasyView, it must be configured with an IP address, this is achieved by using the Device Manager.

Device Manager is located in **Start/All Programs/MKS RGA Applications/Tools/Device Manager**.

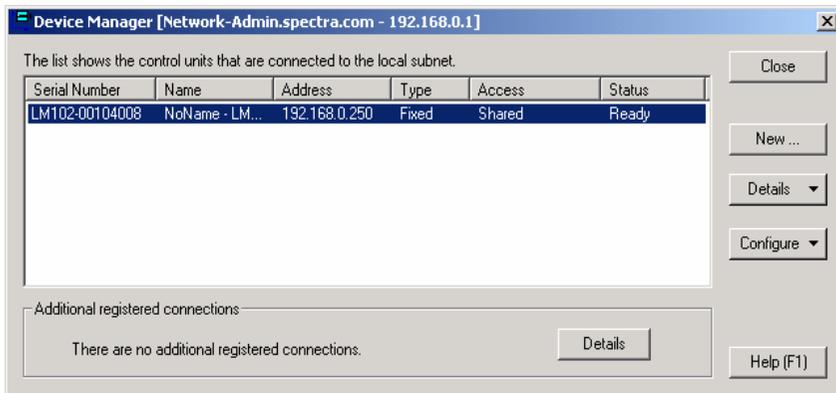

Information pane

"Serial Number" – the serial number of the instrument

"Name" – the friendly name of the instrument

"Address" – the IP address of the instrument

"Type" – whether the IP address is fixed (manually assigned), or DHCP (assigned by your DHCP server) – **this is not supported**

"Access" – see "Access Levels" on page 37.

"Status" – the current status of the instrument

Controls

<Close> - closes the Device Manager dialog

<New> - specify an instrument by its serial number

<Details> - "Auto-discovery log" - displays information on all found sensors

"Refresh" – update the list of discovered sensors

<Configure> - "Browse to the instrument" – displays the homepage of the sensor


"Change IP address" – change to the IP address and subnet

"Access" – the access level of the sensor


<Details> - display a list of past connections that are no longer active

These functions are described in more detail beginning on page 33.

When Device Manager is started, the program searches for connected instruments and displays information on any found, which exist on the same IP scope as the host PC. The host PC's IP address and network name is also displayed in the title bar of the dialog.

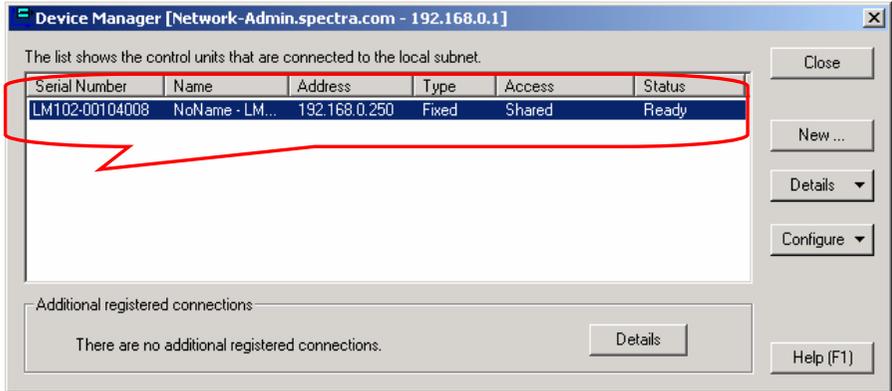

If no instrument information is displayed, it may be that your network uses a scope outside the 192.168.0.0-255 range. In this case, you can use the <New> button and search for the instrument by its serial number.

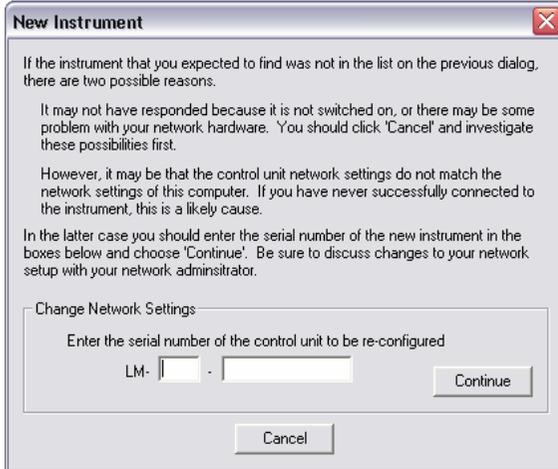


Enter the serial number of the instrument you wish to configure in the text box and click the <Continue> button.

A dialog is displayed warning of the consequences of IP changes to an instrument.

Once your instruments have been installed and are in use on your network, you should observe the warning information displayed, but during this initial configuration, click the <Yes> button to continue.

IP Configuration dialog

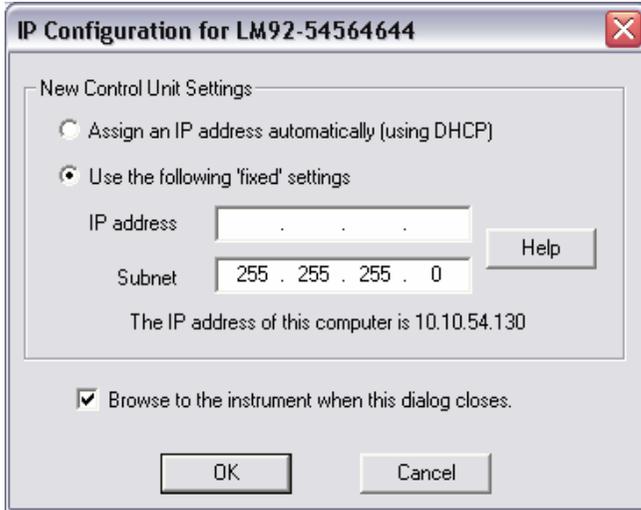

Options

“Assign an IP address automatically using DHCP”
This feature is not currently supported.

“Use the following fixed settings”
The IP address you will assign to the instrument.


“Browse to the instrument when this dialog closes”
Once the address change has completed, your browser is started and you are taken to the instrument's home page. See Section 15 for more information on the web functions of the e-Vision.

“Help”
Launch the help program on IP addressing and networking.

Enter the new IP address of the instrument and click the <OK> button to commit the changes.

The instrument information is now displayed, if no instrument has been found, check that power has been applied and the network cable is attached.

For further troubleshooting advice, see Section 17 - Troubleshooting

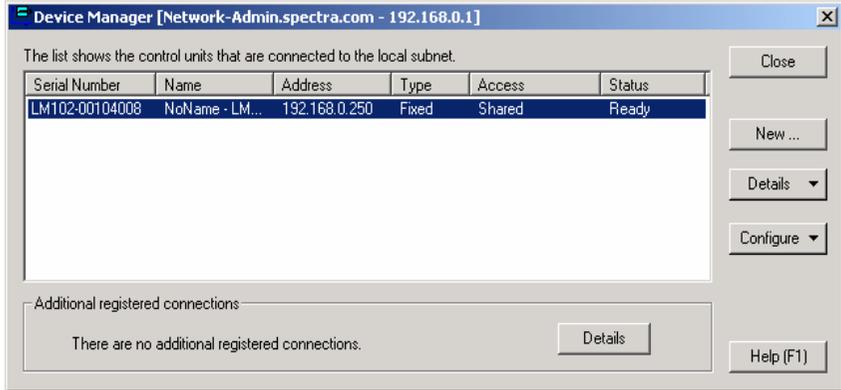

Device Manager Controls

<Details>

“View auto-discovery log” – displays the action history of Device Manager when searching for instruments.



“Refresh” – refresh the current list of found instruments.

<Configure>

"Browse to this instrument" – starts your web browser and takes you to the home page of the instrument.

For more information on the web applet, see Section 16 – Java Web Applet

"Change the IP address" – enter a new IP address for the instrument.

"Additional registered connections"

Click on the <Details> button to display a list of previously registered connections to instrument which are no longer available.

Once all the heads have been configured from Device Manager the installation is complete.

"Access" – set the access level of the instrument, see next page for explanation.

Once the access levels have been configured, you can now start using the EasyView program, see Section 6 – Using EasyView

Access Levels

The “access level” of the instrument determines whether the instrument is available to other instances of EasyView on other computers, or used exclusively by its host PC.

These options are explained below:



EasyView will not auto-connect to the instrument.
Choose this option if the instrument “belongs” to another user running EasyView on another computer.




EasyView will auto-connect and will allow other users to use the instrument when not in use by this PC.




EasyView will auto-connect and will not allow access to other users whether the instrument is in use by its host or not.

The access level of a particular instrument is reflected in the EasyView Loader screen as shown.

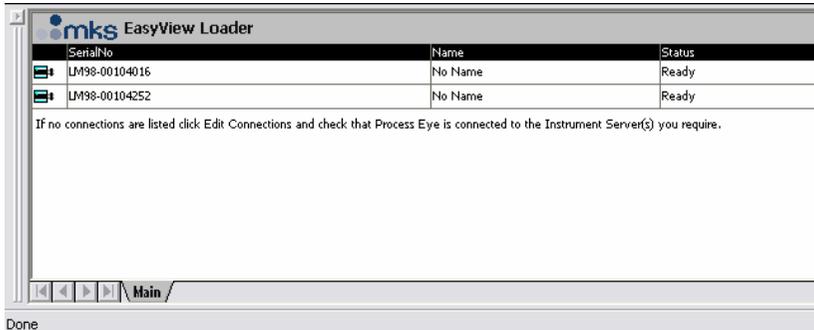

The view above shows two "shared" instruments, EasyView has not made a connection to these instruments, so both are available. You can see that both are "ready" to be used by Easyview. To begin using one of them, highlight it and click on the <Connect> button

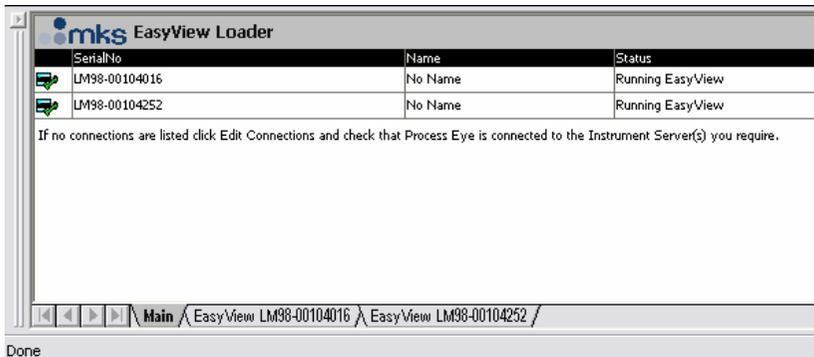

The view above shows two "exclusive" instruments, EasyView has made the connection to both these instances when the program was started. Notice the green tick on both the instrument icons and "running EasyView" is indicated under "Status".

Notice also that both instruments have a "view tab".

6. Using EasyView

Overview


All the standard view modes are available including Analog Scan, Barchart, Peak Jump and Leak Check. These modes are fully interactive allowing changes to the scanning parameters to be made in real-time.

There is also the ability to perform a Disk Store, where the scan data is saved to disk for viewing at a later date using the Recall program.


The following pages describe the main areas of interest to get you started using EasyView. The later sections, describe each area in greater detail as well as explaining the Diagnostics and Calibration modes available.

6.1 Starting EasyView


To start the main EasyView program, click on **Start/Programs/MKS RGA Applications/EasyView**.

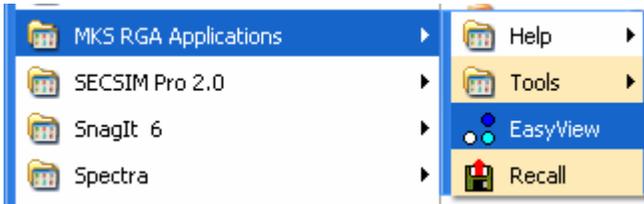

Or by clicking on the desktop shortcut
The CD is not required to run EasyView.

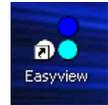

The EasyView information page now loads giving information on the currently connected head.

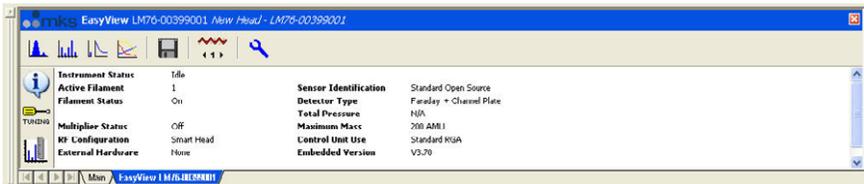

6.2 Toolbars

Each mode of operation can be accessed through the tool bar.



Use the following buttons to start or stop a scan mode, or to access the other functions of the program. For detailed information on each function, see the relevant section shown in brackets.



Analog scan mode. (7.3)



Barchart scan mode. (7.4)



Leak Check scan mode. (7.5)



Peak Jump scan mode. (7.6)



Disk Store. (7.7)



Filament Select.

Use this button to toggle between the filaments.



Filament Status.

Red – Filament ON, Blue – Filament OFF.



Run Diagnostics.

Performs a diagnostics check on the RGA head.



Show Information Page.

This button switches from the Status display to the Information display.



Tuning Mode.

Enter the instrument tuning mode.




Calibration Mode.




Peak Jump Settings.
Only visible when in Peak Jump mode.

6.3 Analog Mode

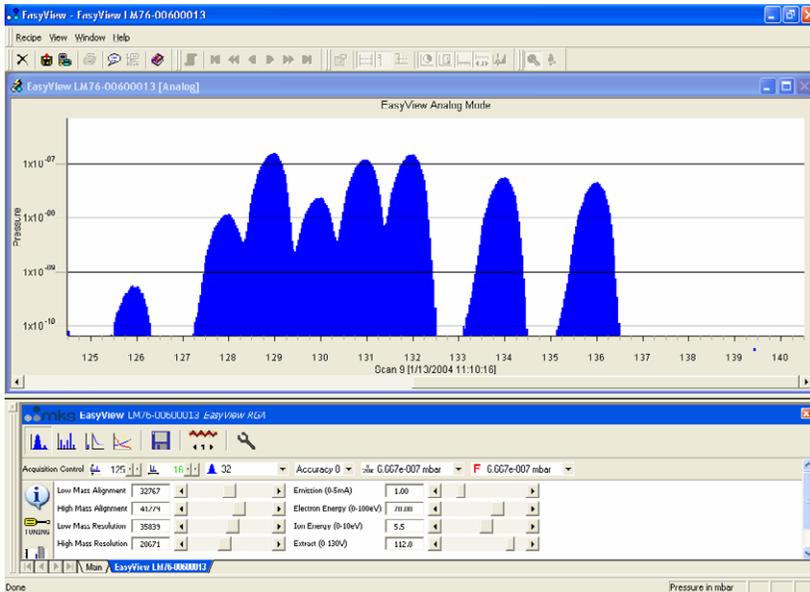

The Analog view is displayed and the interactive analog control toolbar will appear in the status window.

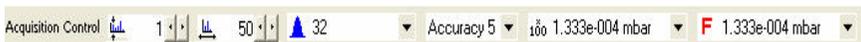

Use this toolbar to change the scan parameters, each of the functions are explained below.



Start Mass control.

To change the start mass of the scan, either enter a number directly into the box, or click on the arrows to increase or decrease the mass number. Note that the first mass cannot be less than 1 and that the first mass plus the mass span, cannot exceed the maximum mass allowed by the control unit.



Mass Span control.

To change the mass span of the scan, either enter a number directly into the box, or click on the arrows to increase or decrease the mass span. Note that the first mass plus the mass span cannot exceed the maximum mass allowed by the control unit.

Accuracy 5 ▼

Accuracy Control.

Scan accuracy determines the amount of averaging carried out for each reading. A high accuracy number provides a clean scan but increases scan time. The accuracy ranges available are from 0 to 8.

▲ 32 ▼

Points per peak control.

The number of measurements each full mass is divided into 8,16 or 32. The higher the number of points-per-peak selected, the better the representation of peak shape. However, scan time and the amount of data generated will increase. If the data is to be imported into Excel for instance, choose a lower value.

100 1.333e-004 mbar ▼

Electronic Gain control.

The electronic gain of the pre-amplifier can be changed by selecting one of the values from the drop down list.

Depending on the type of instrument, the allowed values are 1,100 and 20000. The higher the gain selected, the smaller the partial pressure that can be measured, this may cause peaks of a higher partial pressure to saturate.


F 1.333e-004 mbar ▼

Detector control.


The detector type used for the scan can be changed.

Depending on the type of instrument, the options are Faraday, Multiplier 1, Multiplier 2 and Multiplier 3. Multiplier 1, 2 and 3 refer to the same physical detector, but using three separate calibration values.

6.4 Barchart Mode

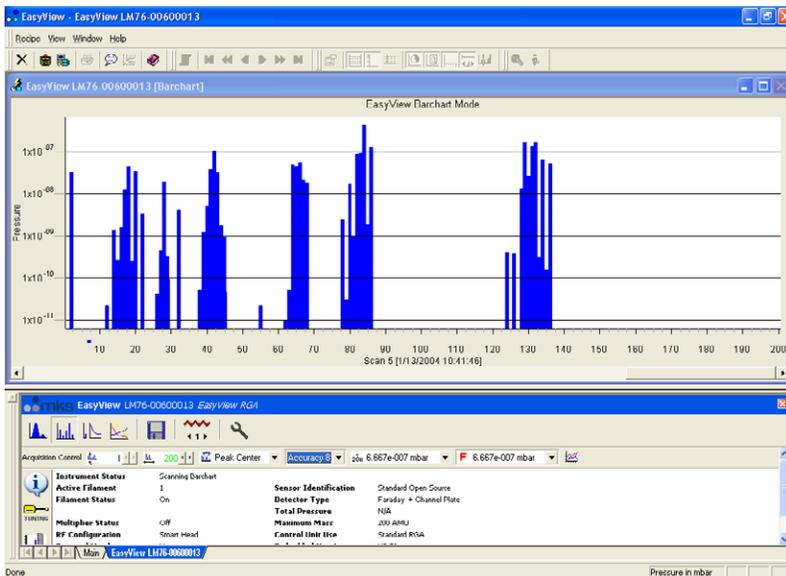

The Barchart view is displayed and the interactive Barchart control toolbar will appear in the status window.

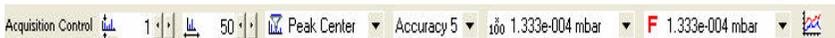

Use this toolbar to change the scan parameters, each of the functions are explained below.



Start Mass control.

To change the start mass of the scan, either enter a number directly into the box, or click on the arrows to increase or decrease the mass number. Note that the first mass cannot be less than 1 and that the first mass plus the mass span, cannot exceed the maximum mass allowed by the control unit.



Mass Span control.

To change the mass span of the scan, either enter a number directly into the box, or click on the arrows to increase or decrease the mass span. Note that the first mass plus the mass span cannot exceed the maximum mass allowed by the control unit.

 Peak Center ▼ Scan Mode control.

The scan mode is the method used to report the peak height.
The options are - Peak Centre, Peak Max or Peak Average.

Accuracy 5 ▼ Accuracy control.

Scan accuracy determines the amount of averaging carried out for each reading.
A high accuracy number provides a clean scan but increases scan time.
The accuracy ranges available are from 0 to 8.

100 1.333e-004 mbar ▼ Electronic Gain control.

The electronic gain of the pre-amplifier can be changed by selecting one of the values from the drop down list.
Depending on the type of instrument, the allowed values are 1,100 and 20000.
The higher the gain selected, the smaller the partial pressure that can be measured, this may cause peaks of a higher partial pressure to saturate.


F 1.333e-004 mbar ▼ Detector control.

The detector type used for the scan can be changed.
Depending on the type of instrument, the options are Faraday, Multiplier 1, Multiplier 2 and Multiplier 3. Multiplier 1, 2 and 3 refer to the same physical detector, but using three separate calibration values.

 Add Trend control.

The add trend button adds a default trend in addition to the Barchart view.
This trend can be edited in the same way as any normal trend view see Section 7 – Changing/adding Views

6.5 Leak Check Mode

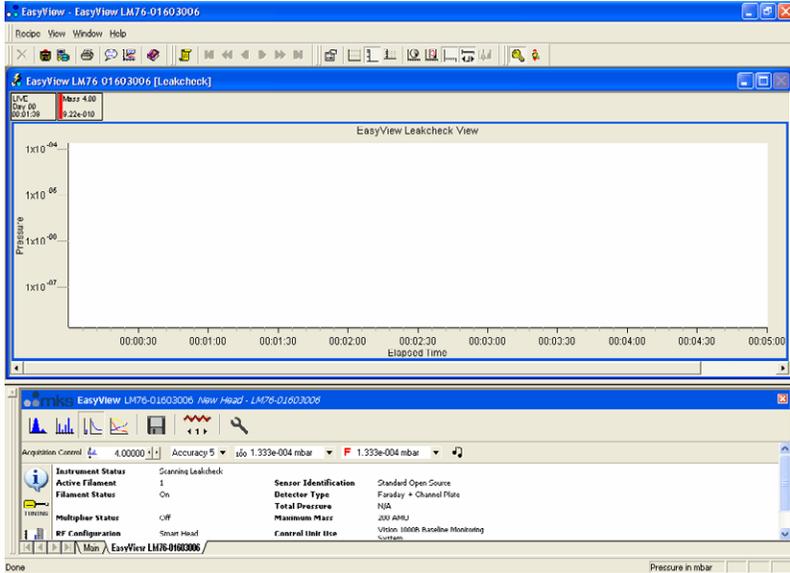

The Leak Check view is loaded and the interactive Leak Check control toolbar appears in the status window.



Use this toolbar to change the scan parameters, each of the functions are explained below.



To change the probe mass of the scan, either enter a number directly into the box, or click on the arrows to increase or decrease the mass number. Note that the first mass cannot be less than 1 and that the first mass plus the mass span, cannot exceed the maximum mass allowed by the control unit.



Scan accuracy determines the sampling amount carried out on each reading. A high accuracy number provides a clean scan but increases scan time. The accuracy ranges available are from 0 to 8.

100 1.333e-004 mbar ▼

Electronic Gain control.

The electronic gain of the pre-amplifier can be changed by selecting one of the values from the drop down list.

Depending on the type of instrument, the allowed values are 1,100 and 20000. The higher the gain selected, the smaller the partial pressure that can be measured, this may cause peaks of a higher partial pressure to saturate.

F 1.333e-004 mbar ▼

Detector control.

The detector type used for the scan can be changed.

Depending on the type of instrument, the options are Faraday, Multiplier 1, Multiplier 2 and Multiplier 3. Multiplier 1, 2 and 3 refer to the same physical detector, but using three separate calibration values.




Audio control. Not support on certain sensor configurations.

Switches the audio tone on or off.

The tone changes in frequency as the measured peak height changes.

This is a useful feature if you are unable to see your monitor while leak checking.

6.6 Peak Jump Mode

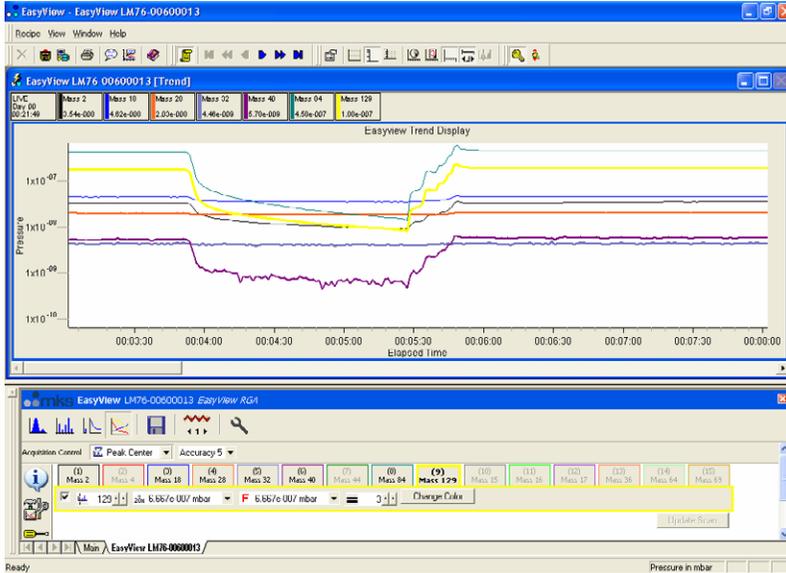

The Peak Jump view is loaded and the interactive Peak Jump control toolbar appears in the status window.



Use this toolbar to change the scan parameters, each of the functions are explained below.




The scan mode is the method used to report the peak height. The options are - Peak Centre, Peak Max or Peak Average.



Scan accuracy determines the sampling amount carried out on each reading. A high accuracy number provides a clean scan but increases scan time. The accuracy ranges available are from 0 to 8.



Use this button to open the Peak Jump settings window shown on the next page.

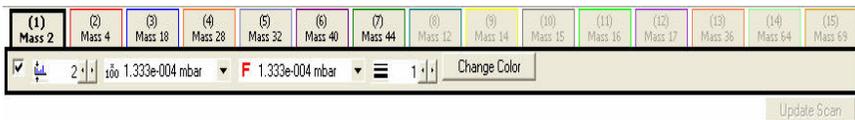

Settings for each of the 15 channels can be edited using this toolbar.




Click on the channels tab to edit the settings for that particular channel.



Check the box to enable the channel.

In the above example only channels 1 to 7 are enabled, channels 8 to 15 are disabled.

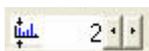

Scan Mass control.

Set the particular channels scan mass.

Note that the mass cannot be less than 1 and cannot exceed the maximum mass allowed by the control unit.



Electronic Gain control.

The electronic gain of the pre-amplifier can be changed by selecting one of the values from the drop down list.

Depending on the type of instrument, the allowed values are 1,100 and 20000. The higher the gain selected, the smaller the partial pressure that can be measured, this may cause peaks of a higher partial pressure to saturate.



Detector control.

The detector type used for the scan can be changed.

Depending on the type of instrument, the options are Faraday, Multiplier 1, Multiplier 2 and Multiplier 3. Multiplier 1, 2 and 3 refer to the same physical detector, but using three separate calibration values.



Line Thickness control.

The thickness of the plotted line can be changed to aid visibility



Colour control.

The colour of the plotted line can also be changed, select a colour then click <OK> to apply it to the channel.




Update Scan

Changes made to the channels properties are not immediately made to the scan. When all changes required have been made click <Update Scan> The new settings will then be applied to the current scan.

6.7 Disk Store




Disk Store is available in any of the EasyView scan modes and can be accessed by clicking the button shown.

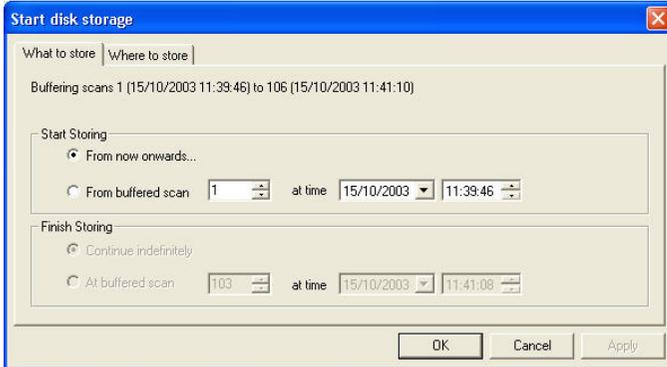

What to store options:

"From now onwards" stores each buffered scan to disk until stopped by the user.
"From buffered scan" allows *start* and *end* scan configuration.

Note: The buffer is reset if any changes to the current scan are made, otherwise the last 500 scans are stored in the buffer.
If disk storage is started at a buffered scan, then it is possible to stop disk storage at a defined buffered scan as well.

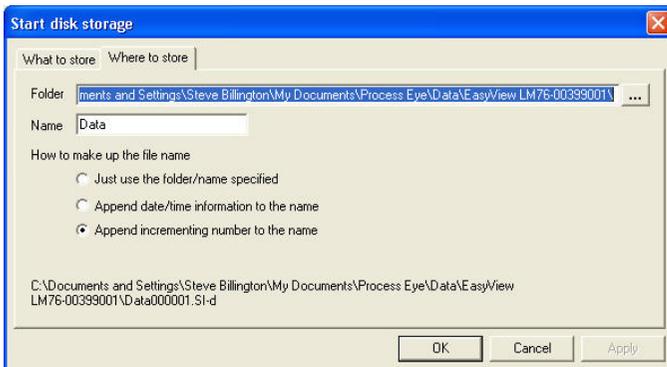

Where to store options:

"Folder" enter the path, or browse to the folder where the scans are to be stored. The default path is:

..\EasyView\Data\EasyView "Head Serial Number"

"Name"

Enter a name for the saved data file, by default this is "data" but it can be anything you choose.

As well as the name, the file can be amended with other information:


"Just use the folder name specified" will create the file "data.SI-d".

"Append date/time information to name" adds date and time information to the file and would create the file "DataYYYYMMDDHHmmSS.SI-d".

Where YYYY is the current year, MM is the current month, DD is the current day, HH is the current hour, mm is the current minute and SS is the current second.

"Append incrementing number to the name" adds a sequential number to the name and would create the following file "Data000001.SI-d", the next would be "Data000002.SI-d" and so on for each file.

While storing to disk, the Disk Store toolbar will be visible in the status window.

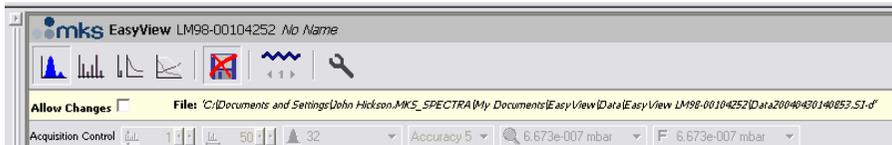

Notice that the disk store icon is no longer available and the path to the disk store file is displayed.

Checking the "Allow Changes" box, pauses the disk store and re-enables the scan toolbar.




You can then change the scan properties, un-check the "Allow Changes" option and continue with the Disk Store without having to run through the "Start Disk Storage" dialog.

Note: The name of the file will still change.

6.8 The Main EasyView screen

The main EasyView screen is where most of your time will probably be spent. The picture shows the main screen in Analog scan mode.

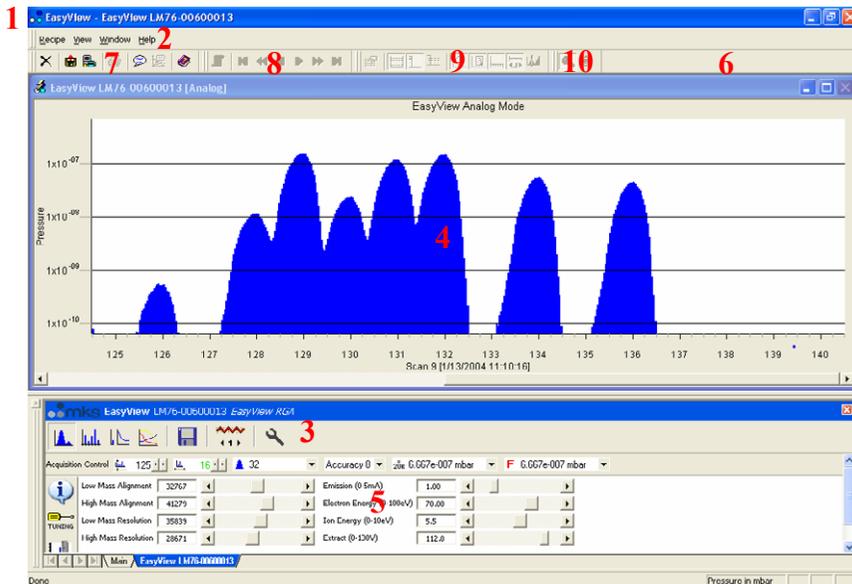

1 Title Bar

The Title Bar shows the application name, EasyView. The Close, Minimize and Maximize button on the right-hand side of the title bar operate in the normal manner.

2 Menu Bar

EasyView is no different to other Windows programs, all the available function are available via the menu bar.

3 EasyView Control Icons

EasyView can be controlled from these icons, use them to select the required mode.

4 Views Windows

The Views windows are where the data acquired in EasyView is displayed. Any number of trend windows many be added to the default views.

5 EasyView Control Window

The EasyView control window contains all the controls to run EasyView. From here select the scan mode, control disk store, set the control unit parameters, calibrate and run diagnostics

6 Toolbars

The Toolbars appear as one row of buttons below the menu bar and immediately above the recipe list and views windows. The twenty-five buttons consist of four separate toolbars. Each of the toolbars may be hidden by selecting <View / Toolbars> from the menu bar and un-checking those you wish to hide. The default setting is visible for all four toolbars.

7 Default Toolbar



The first seven buttons make up the Default Toolbar.



The "Close" recipe button is used to halt a recipe that is currently running. The Close Recipe button will not be available unless a recipe is running and the Status window has focus (click in the status window to give it focus). Also, the recipe needs to have been configured to include the ability to end the recipe prematurely.



The "Run Recall" button will run the recall program to allow review of stored data



The "Run Software Setup" button will launch the software setup wizard to allow changes to the installation



The Print button will print the data from the active data window.



The "Add Annotation" button can be used to add annotation to a trend at a particular time.



The add annotation at selected time button is used to mark a place on a trend



Click on the "Help" button to open the EasyView help file. Use EasyView Help just as you would any other Windows help document.

8 Scan Navigation Toolbar



These seven buttons control the viewing of live and stored data.



When the Historic Data button is not depressed, the active continuous mass scan view is displaying the live data currently being acquired. While the button is depressed, the continuous scan window will display historic data. New data is still being acquired in this historic mode, you will see the total number of scans figure increasing while the current scan number remains the same.

The six navigation buttons will only become available when historic data is being viewed, e.g. The Live Data button is not depressed.



Click on the "First Scan" button to view the oldest scan stored in memory.



Click on the "Jump Back" button to view the scan 10 scans previous to the current one being view.



The "Previous Scan" button allows you to view the last scan.



The "Next Scan" button allows you to view the next scan.



Click on the "Jump Forward" button to view the scan 10 scans subsequent to the current one being view.



Click on the "Last Scan" button to view the newest scan acquired in the current recipe run.

9 Chart Display Toolbar



These nine buttons that make up the Chart Display Toolbar control certain aspects of the active recipe windows.



Click on the "Properties" button to display the settings for the currently active recipe view.



Click on the "Grid" button to display a Y-axis grid in the currently active view window. Default is disabled.



Click on the "Log/Lin" button to switch between a linear and logarithmic Y-axis. Default is linear.



Click on the "Y-axis Minor Ticks" button to enable the display of minor ticks on the Y-axis. Default is major ticks.



Click on the "Time" button to switch between real and elapsed time on the X-axis from the start of the recipe. Default is elapsed time.



Click on the "Days" button to display the date and time on the X-axis. Default is time only.



Click on the "X-axis Minor Ticks" button to enable the display of minor ticks on the X-axis. Default is major ticks.



Click on the "Add Scroll Bar" button to add a scroll bar to the X-axis. Default is no scroll bar.



Click the "Add Cursor" button to add a measurement cursor to the active Bar chart display. The cursor can be moved to the required mass and will display the mass number and partial pressure. As many cursors as are required can be added.

10 Trend Display Toolbar



These two buttons constitute the Trend Display Toolbar



The "Key Display" button controls the display of the key table for the active trend views. The button will not be available if no trend view is available.




The "Key Mode" button changes the key table view between vertical and horizontal views.

Customizing Toolbars

You can customize the toolbars as required, you can choose which Toolbars are visible, or you can create your own toolbars with your own choice of buttons.


From the view pull down menu, select <Toolbars>.

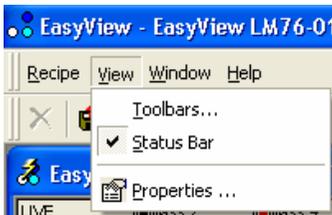

This will bring up the toolbar dialog. Check the box for the toolbars you wish to have displayed. Checking the <Cool Look> check box will change the appearance of the toolbars.




To create a new toolbar, click on the <New> button.

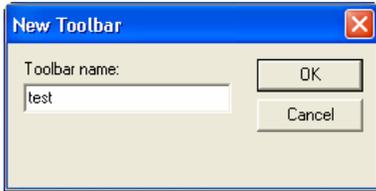

Give the new Toolbar a name and click <OK>.

This will create a blank Toolbar that you can then customize. From the toolbar dialog box click on the <Customize> button to bring up the customize dialog.

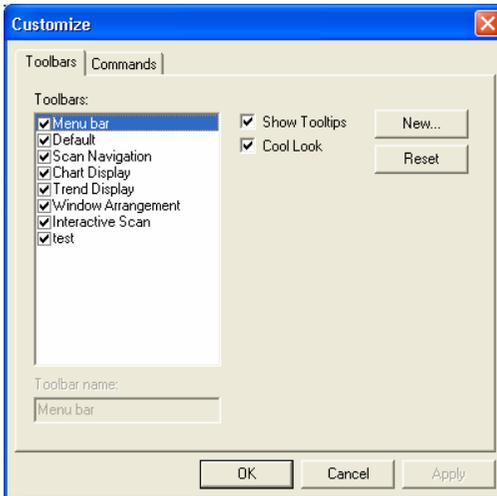

Click on the <Command> tab to see the buttons currently assigned to the new Toolbar you have created.

Adding Annotations

Time stamped annotations can be added to a trend in two ways:

The first is to click on the "Add Annotation" button which will cause the "Annotations" dialog to appear.




The "Annotations" dialog box has a title bar with a close button. It contains a "Time of annotation:" field with a date dropdown set to "28/11/2002" and a time dropdown set to "14:48:02". Below this is a "Note:" text area with a scroll bar. At the bottom are "OK" and "Cancel" buttons.

Here you can specify a time for the annotation to appear on the trend, add the text you wish to appear in the "Notes" field.

The second way to add an annotation is to click on the "Add annotation at selected time" button.



The screen cursor will change to a pen, select the point on the trend that you want to add the annotation and click.

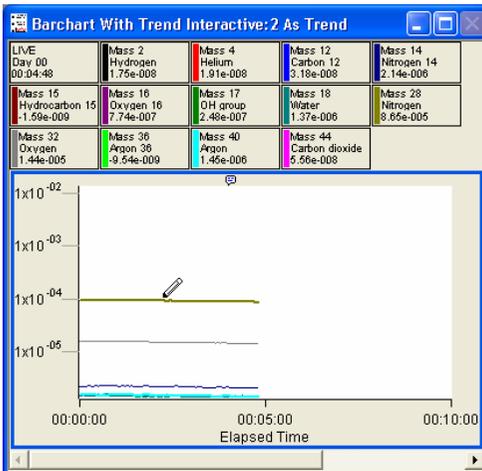

The annotations dialog appears again, allowing you to add text to the annotation, with the date and time determined by the point at which the annotation was added to the trend.



Each annotation added is marked by a symbol along the top of the trend, moving the cursor over this symbol causes the entered text to be displayed as a tool tip. You may add as many annotations as you wish, these will also be visible when using Recall.

7. Changing/Adding Views in EasyView


The RGA data gathered by the instrument can be displayed on screen in a number of different views.

To modify any view manually, highlight the view and either right-click, or click on the properties button on the chart display setup toolbar.




View types


“Swept Measurement” view, this is a display of partial pressure versus mass. Bar chart shown below, where there is a single bar shown per mass.

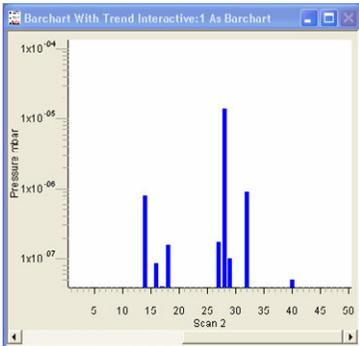

Another form of the swept measurement view is the Analog view, where there are a number of bars per mass, giving a pseudo-analog view.

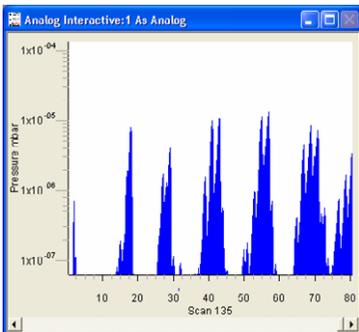


"Trend" view where a variable versus time can be plotted.

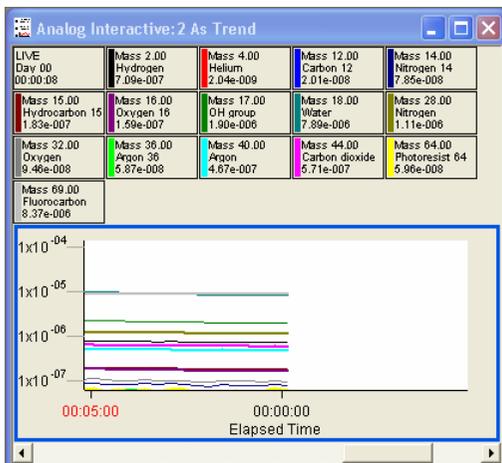

This type of view is more flexible as any user variables can be plotted.

7.1 Swept Measurement Properties Settings

The following section outlines the various changes that can be made to views using the "Properties for Continuous Mass Scan View" dialog. We will look at each function in turn.

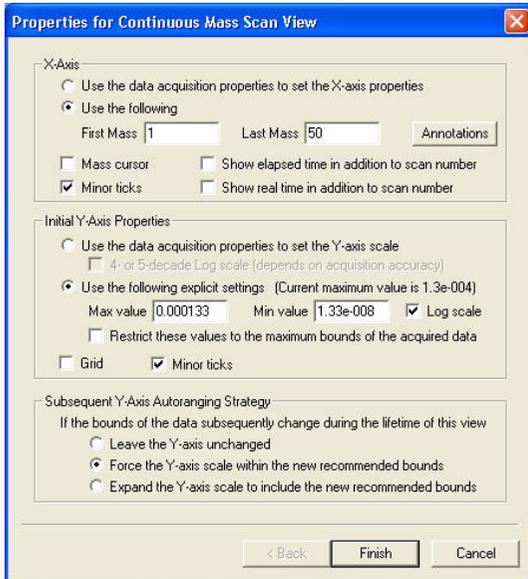

X-Axis

Use the data acquisition properties to set the X-axis properties

Creates a view with the X-axis properties governed by the measurement on which the view has been based.

For example, if the view was based on a Bar chart scan, measuring from mass 5 up to mass 45 then this view will show mass 5 to mass 45 along the X-Axis.

Use the following

Bases the X-axis range on the values entered in the text boxes, with the following limitation:

Although you can choose not to display masses that are being measured, you cannot choose to display masses that are not being measured.

For example, you create a Bar chart measurement scanning from mass 5 to 45, you can base a view displaying from mass 10 to 40, but not from mass 5 to 50. If this is attempted, the view will revert to displaying the measurement's range.

Minor Ticks

Displays minor ticks on the X-axis.

Show elapsed time in addition to scan number

Displays the elapsed time from the start of the recipe.


Show real time in addition to scan number

Displays the real time, governed by the PC clock.

Note: Both the above options cannot be enabled at the same time.

To edit the Annotations icon click on <Annotations>

Annotations dialog

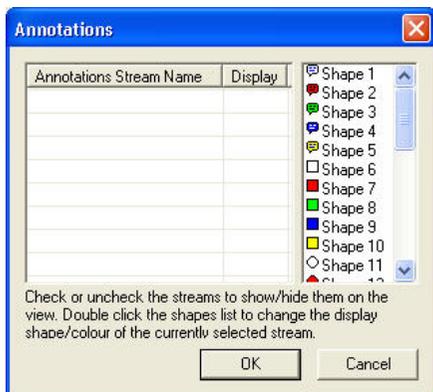

Any annotation streams defined in the recipe will be listed, check the display box to show the annotation on the view. To change the icon displayed on screen, select one of the others from the list.

Click <OK> to return to the properties page

Initial Y-axis Properties

In a swept measurement, the Y-axis is always shown as pressure, the units of which are determined in the EasyView preferences, Section 13 - Preferences.

Use data acquisition properties to set Y-axis

Creates a view with the Y-axis properties governed by the measurement on which the view is based.

4- Or 5- decade log scale

Changes the Y-axis from a linear scale with its origin at zero, to a logarithmic scale, with a minimum on which is dependant on the accuracy of the measurement. The higher the accuracy, the lower the minimum pressure will be displayed.

Use the following explicit settings

Allows manual entry of the maximum and minimum displayed pressures. The display can be either linear or logarithmic determined by the check box.

Restrict these values to the maximum bounds of the acquired data

Causes the Y-axis maximum pressure setting, to be governed by the maximum pressure obtained from the measurement associated with the view.

Note: You cannot display zero or negative values in logarithmic mode.

Grid

Displays a horizontal grid, aligned with the ticks of the Y-axis.

Minor Ticks

Displays minor ticks on the Y-axis.

Subsequent Y-Axis Auto-Ranging Strategy

If a view's Y-axis limits are based on an "interactive" measurement, these limits can be changed during the lifetime of the view.

Three options determine how the axis will react to changes in the measurement such as electronic gain range, or detector changes for example.

Leave the Y-Axis unchanged

The Y-Axis settings remain unchanged when the measurement changes. The settings will be based on the measurement when the view is loaded.

Force the Y-Axis scale within the new recommended bounds

Changes the Y-Axis settings to match the new measurement settings.

Expand the Y-Axis scale to include the new recommended bounds

Change either the maximum or the minimum Y-axis settings to incorporate the new settings.

For example:

If you increase the electronic gain, the minimum displayed pressure would be changed to match the smaller pressure that could now be measured. The maximum pressure displayed would remain the same. If the electronic gain was then reset to its previous value, the Y-Axis would not change.

Once the changes to the view settings are complete, the view can be saved. Click on the <Next> button to bring up the "Save View" dialog box.

7.2 Trend View Properties Settings

Trend views

A trend view can be associated with any measurement. Along with the RGA data, it is also possible to display signals from the analog / digital inputs and other user variables. The only limitation is that the X-axis is always time.

Adding a Trend view

To add a trend view simply click on the new trend button on the EasyView control page.




Alternatively select <New Trend Window> from the <Window> pull down menu



The first dialog displayed is the "Trends Display Pane and Channel Properties". By default, 13 of the 15 masses are enabled, but you can enable and disable channels by checking the box next to the channel.

Clicking on the "Use standard view" will set the channels to be the same as those in the default trend.

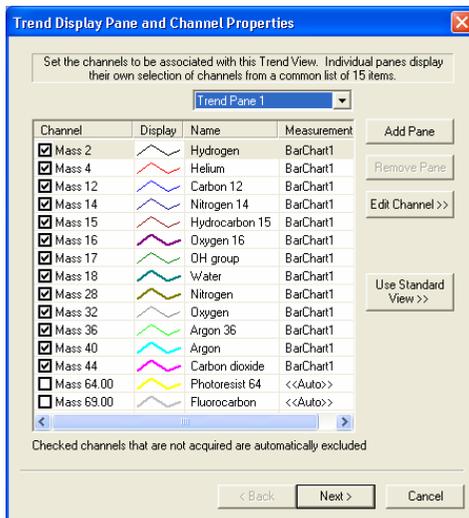

The list of common gases can be edited to suit your needs, see Section 12 – Preferences for details.

Editing a Trend channel

You can edit a channel by either double-clicking it, or by highlighting the channel and then clicking the “Edit channel button”.

This displays the “Channel Properties” dialog, which allows set-up of each channel independently.

The Channel Properties options are explained below.

Channel Properties

Entry for channel 1
Type: Mass Value: 2.00

Optionally enter a measurement name, or leave blank for auto-selection when this view is used

Channel Name or Alias: Hydrogen

Color: [Black] Line thickness: [Slider] Draw Stepped:

For mass channels you may double click an item from the list

Mass	Name
1	Atomic hydrogen
2	Hydrogen
4	Helium
12	Carbon
14	Nitrogen(2)
14	CH2 group
14	Atomic nitrogen
15	CH3 group
15	Methane(2)
16	Methane
16	Atomic oxygen

OK Cancel

Type

“Mass” refers to one of the partial pressure measurements.

Type: Mass

Editing a mass channel

If "Mass" is selected as the channel type there are a number of ways of editing the channel.

One way is to select one of the pre-defined channels in the selection table.

Mass	Name	
1	Atomic hydrogen	
2	Hydrogen	
4	Helium	
12	Carbon	
14	Nitrogen(2)	
14	CH2 group	
14	Atomic nitrogen	
15	CH3 group	
15	Methane(2)	
16	Methane	
16	Atomic oxygen	

Clicking on one of the lines will set the mass value and channel name.

Alternatively, you can type in the value of the mass you wish to display in the "Value" text box.

Then type the name that you wish to appear on the trend key in the "Channel name or alias" text box.

If you leave this box blank, the channel name will be displayed as "Mass ??.??" where "??.??" is the value entered in the "Value" text box.

The measurement name is the name of the measurement that you want to use to provide the partial pressure data. If the box is left blank, the measurement selection will be "Auto".

This does not mean that the best data will be selected, only that the software will look for a mass with the value entered for each measurement in the measurement list. The first measurement that contains this mass value will be the one that is used.

For example:

You have two Bar chart measurements, Bar chart 1, that is a mass 1 to 50 Faraday measurement and Bar chart 2, which is a mass 1 to 100 multiplier measurement.

If the channel has a mass value of 4.00, the data would be taken from Barchart1 because this is the first measurement that contains a mass with a value of 4.00.

This will be true, even if the peak height on mass 4 is so small that it is lost in the noise level and mass 4 in multiplier is a good stable peak.

If the value is mass 84 the data would be taken from Barchart2, even if the peak is so large it is off-scale. This is because Barchart2 is the first measurement that contains mass 84.

If the measurement name entered is not a valid measurement name in the current recipe, then the measurement selection defaults to auto.

Editing the channel display

To change the colour of a trend line, click on the "Colour" list box.



This will bring up a standard Windows colour selection dialog, where you can choose a new colour.

To change the thickness of the displayed line, move the slider on the "Line Thickness" control.



Moving the slider to the right increases the thickness.
Moving the slider to the left decreases the thickness.

The "Draw Stepped" check box can be used when displaying digital information on a trend.




Instead of each point on the trend being joined by a straight line, the points are joined only by horizontal and vertical lines.

The data is forced to be a value of either zero or one. This feature is particularly useful for displaying digital channels on a trend.

Channel	Display	Name	Measurement
<input checked="" type="checkbox"/> Mass 14.00		Atomic nitrogen	<<Auto>>


Adding a Trend pane

Although you cannot define multiple Y-Axis scales to allow display of different data on the same trend, you can create multiple trend panes, which allow you to display a number of different Y-Axis on the same trend.

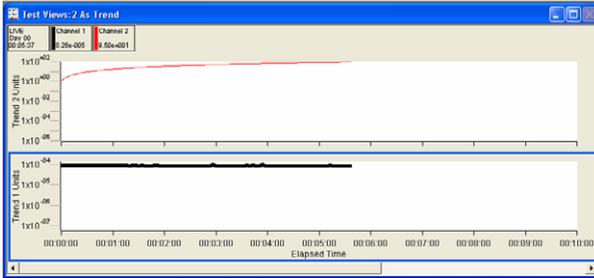


This is a 2-pane trend with channel 1 displayed on trend pane 1, with its own measurement units, and channel 2 displayed on trend pane 2 with different measurement units.

On the "Trend display pane and channel properties" dialog, click on the <Add pane> button to add a new pane.

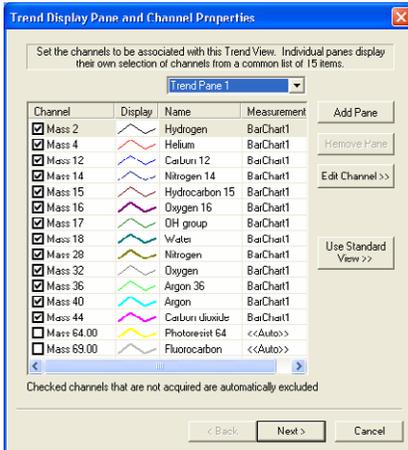

The settings for each pane can be modified. Select the pane that you wish to edit from the Trend Pane Pull-down list box.

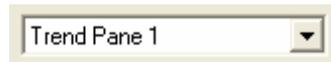

The two-pane view at the start of this section was created in the following way.

Trend Pane 1

Channel	Display	Name	Measurement
<input checked="" type="checkbox"/> User defined		Channel 1	UserMeasure.
<input type="checkbox"/> User defined		Channel 2	UserMeasure.

Channel 1 was enabled on trend pane 1


Trend Pane 2

Channel	Display	Name	Measurement
<input type="checkbox"/> User defined		Channel 1	UserMeasure.
<input checked="" type="checkbox"/> User defined		Channel 2	UserMeasure.


Channel 2 was enabled on trend pane 2

Configuring the Trend's Y-Axis

Once the channels have been set, click on the <Next> button causing the "Properties for Trend View Pane Y-axis" dialog to appear.

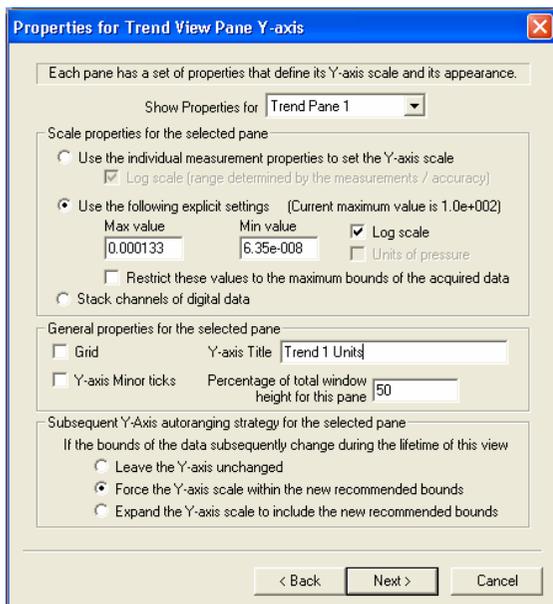

Each trend pane can have a different Y-Axis setup. To set up each pane, select the pane you want to edit from the pull-down list box.

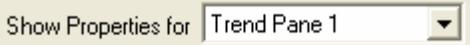

Y-Axis Scale Properties

The scale can be set in a number of ways.



By selecting the above option, the scale will be defined by the measurements that the displayed channels use.

E.g. full scale on the axis will be equal to the full-scale pressure of the measurement.

Checking the “Use the following explicit settings” option allows the maximum and minimum limits on the Y-axis to be defined manually.



Enter the maximum and minimum values you wish to display in the text fields.

Max value
0.000133

Min value
6.35e-008

To display a Log scale, check the box.

Log scale

Note: The software cannot display negative numbers or zero when set to Log scale.

If the “Units of pressure” check box is enabled, the axis will change scale if the pressure units are changed in the preferences.

Units of pressure


Restrict these values to the maximum bounds of the acquired data

By checking the above option, the scale is limited by the maximum limits of the measurements being used.

For example, if the only data to be displayed is mass channels from Barchart1, the limit of the maximum value you can have, will be the full-scale pressure on Barchart1.

With this option selected, the Y-axis does not have a scale as such. All enabled channels are given a value of either zero or one and are then plotted against their channel name.

Stack channels of digital data

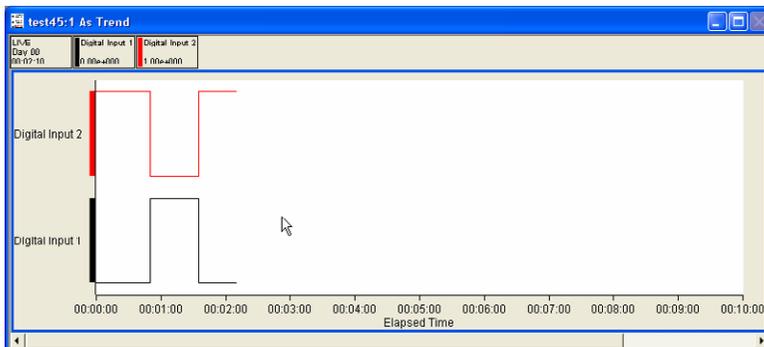

This option is only useful for displaying digital data. In the above example, there are two digital input channels with both channels configured to “draw stepped”.

Y-Axis General Properties

Enabling the grid check box will draw a horizontal grid on the trend.

 Grid

Enabling the minor ticks check box will display minor ticks. If grid is enabled, grid lines for the minor ticks are also displayed.

 Y-axis Minor ticks

Enter a label for the Y-axis in this text box.

Y-axis Title

If there is only one trend pane on the trend this box is set to 100% and unavailable.

Percentage of total window height for this pane

If there is more than one trend pane, you can set the proportion of the trend window that the particular pane will occupy.

Percentage of total window height for this pane

Y-Axis Auto-range Properties

- Leave the Y-axis unchanged
- Force the Y-axis scale within the new recommended bounds
- Expand the Y-axis scale to include the new recommended bounds

If the view uses an interactive measurement to set the Y-Axis limits, then those limits can change during the lifetime of the view. Changing the electronic gain, or type of detector for example.

You can select one of three options to control how the view will react to a change in the measurement:

Leave the Y-Axis unchanged

Leaves the Y-Axis settings unchanged when the measurement changes. The settings will be based on the measurement settings when the view is loaded.

Force the Y-Axis scale within the new recommended bounds

Changes the X-Axis settings to match the new measurements settings.


Expand the Y-Axis scale to include the new recommended bounds

Changes either the maximum or minimum Y-axis setting, to incorporate the new settings.

For example, if you increase the electronic gain, the minimum displayed pressure would be changed to match the smaller pressure that could now be measured. The maximum pressure displayed would remain the same. If the electronic gain was reset to its previous value the Y-Axis would not change.

Configuring the trend time axis

Click <Next> on the Trend properties dialog to bring up the "Time axis and legend properties" dialog.

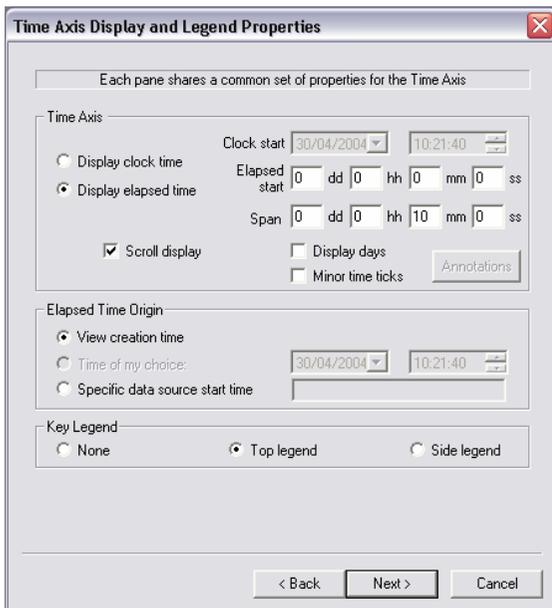

Time axis properties

This option will display the time axis using real time as calculated by the PC clock.

Display clock time

This option displays the time axis as an elapsed time. The origin of the elapsed time will depend on how the time origin is defined (See the next section on configuring the elapsed time origin).

Display elapsed time

Offset from the elapsed time origin that the time axis will start to display.

Elapsed start dd hh mm ss

Amount of time that will be displayed on screen, in this case the last 10 minutes of data will be displayed.

Span dd hh mm ss

Scrolls the Trend to keep the latest data displayed.

If unchecked, the trend will only display data for the first x minutes. Where X is the number of minutes set in the span.

Scroll display


Displays the number of days, as well as the time the Trend has run.

Display days

Displays minor ticks on the time axis.

Minor time ticks

To edit the Annotations icon click on <Annotations>

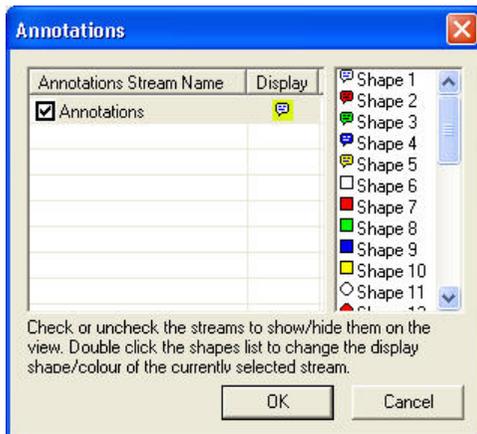

Any annotation streams defined in the recipe will be listed, check the display box to show the annotation on the view. To change the icon, select one of the icons from the list.

Click <OK> to return to the properties page

Elapsed time origin properties

Displays the creation time of the view.

View creation time

This option is only available when editing a live trend.

Time of my choice:

Sets the elapsed start time, to the same time that the measurement was created. Type the name of the measurement in the text box.

Specific data source start time:

Key Legend properties

None – No legend is displayed

Top Legend – Display the legend at the top

Side Legend – Display the legend at the side

Editing a live Trend

Trend views can be changed by using the following toolbar buttons, these only affect the currently active view.

Alternatively, right-clicking the Trend will show the "View Properties" dialog.



Click the Y-axis "minor ticks" button to turn minor ticks on or off for the Y-axis.



Click the Y-axis "grid" button to turn the grid on and off for the Y-axis.



Click the Y-axis "log/lin" button to change between log and linear mode for the Y-axis.



Click the "display days" button to switch on and off the "display days" function.



Click the "true/elapsed time" button to switch between displaying the "real time" and the "elapsed time".



Click the "X-axis minor ticks" button to turn minor ticks on and off for the X-axis.



Click the "X-axis scroll bars" button to turn scroll bars on and off for the X-axis.




Click the "legend key" button to make the legend key visible or invisible.



Click the "table view" button to move the legend key either to the left of the trend or above the trend.


Adding annotation to a Trend

To add annotation, click the <Annotation> button, this will bring up the "Annotations" dialog box.

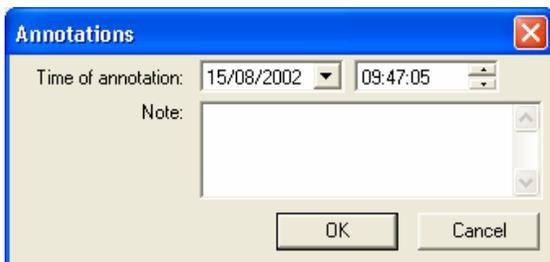

Here you can enter text into the "Note" text box, when your annotation is complete, press the <OK> button. By default, the time of annotation will be the current time, although you can select a new time, by using the "time" and "date" list boxes.

To add an annotation to a trend, click the <Trend Annotation> button, this will change the cursor to a pen.

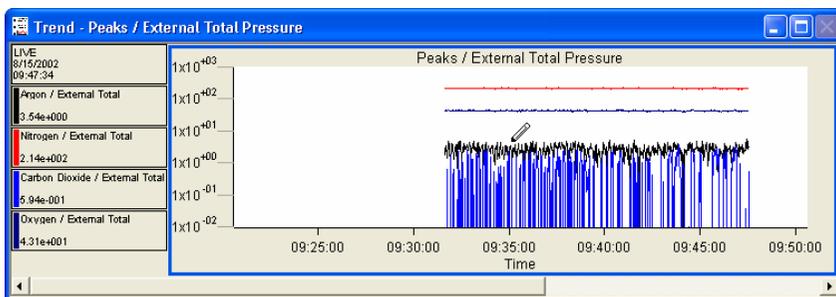

You can move the cursor along the trend line, until the event that you want to annotate, has been selected. Clicking on the trend at this point will bring up the annotation dialog, but now the time will be set to the time at that position on the trend.


Currently, the only way to view these annotations is to recall a saved file with annotations, and then save this file as a text file.

The annotations cannot be viewed in any of the EasyView programs.


Once added, annotations cannot be removed as they become part of the measurement data.

View Global Properties

You can set the global properties of the views from the “Graph Settings” under “Preferences”.

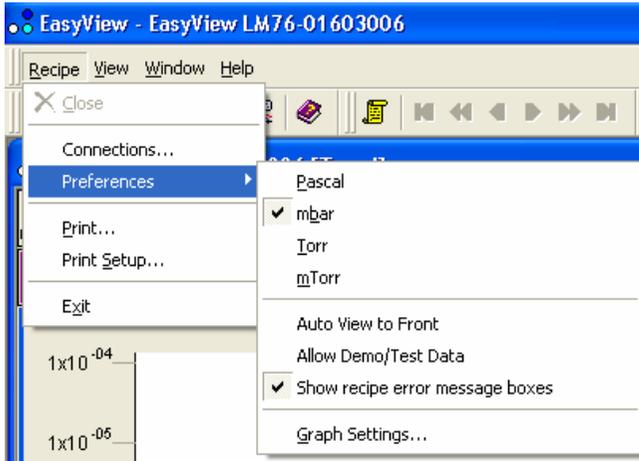

The “Graph Preferences” dialog will appear.

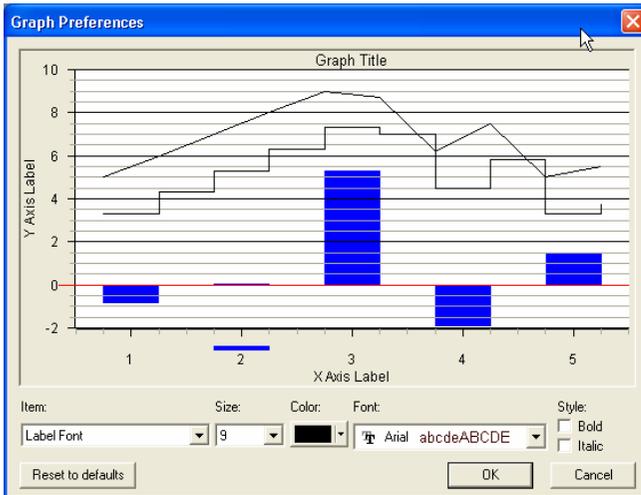


Here you can manipulate the various options used in the graphs, size, colour, font and the style of each item.

The dialog changes to reflect any changes that are made.


Clicking on the <Reset to Defaults> button will return all the settings, to their factory defaults.

The following diagram shows all the options that can be set. If the properties cannot be changed for a selected item, they will remain "greyed-out".

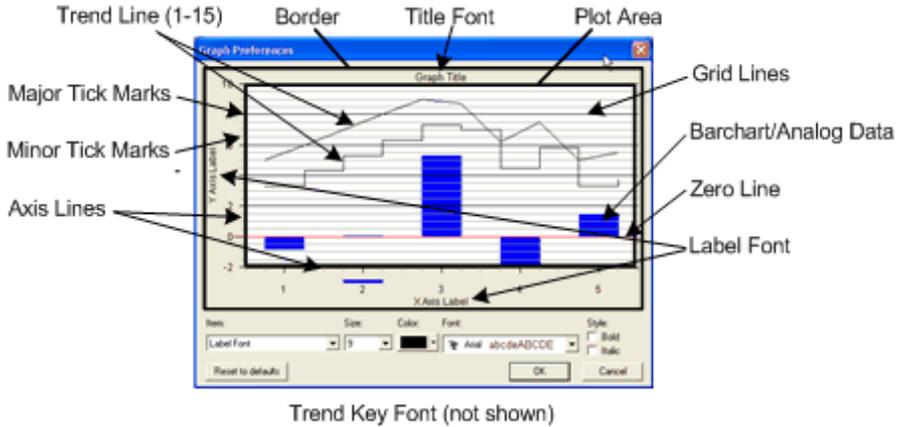

8. Other Menus

View Pull-Down Menu

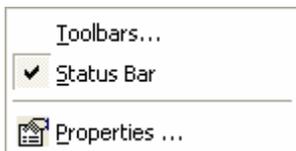

This menu can be used to set up the appearance of EasyView.

Toolbars

Selecting "Toolbars" from the menu, will bring up the Toolbars edit dialog.



Status Bar

The Status Bar appears at the bottom of the screen.



To make the Status Bar visible/invisible, click on the "Status Bar" item in the pull-down menu. If a check is visible, then the Status Bar will be visible.

Properties


The Properties item will only be active if a mode is running with a view loaded.

Click on the properties button to bring up the view properties of the active view. If the view is a trend, you will get the trend properties dialog

If the view is a swept measurement view then you will get the swept measurement view properties dialog.

Window pull-down menu

The "Window" pull-down menu is where you can add a new trend view or arrange the view windows you already have.

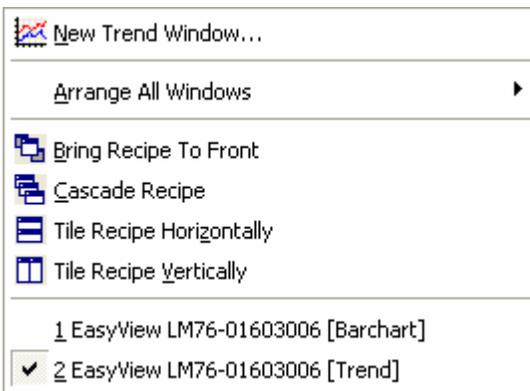

New Trend Window

Clicking on new trend window will bring up the trend properties window for a new view.

Arrange All Windows

This item will bring up an additional list.



From here, you can arrange the windows in the usual way. In this menu however, all the currently open views from all the currently running measurements will be arranged.


The other arrangement methods in this pull down menu, apply only to the active recipe.

Bring Recipe to Front

Selecting Bring Recipe to front will bring the views for the current active measurement to the front


Cascade Recipe

Selecting cascade recipe will bring the views for the currently active measurement to the front and arrange them in a cascade fashion.

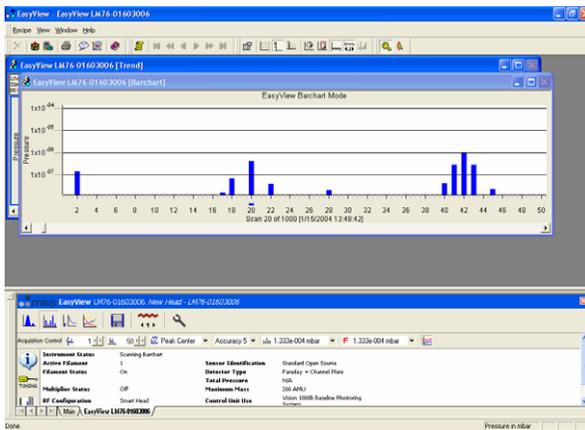


Tile Recipe Horizontally

Selecting tile recipe horizontally will bring the views for the current active measurements to the front and tile them horizontally.

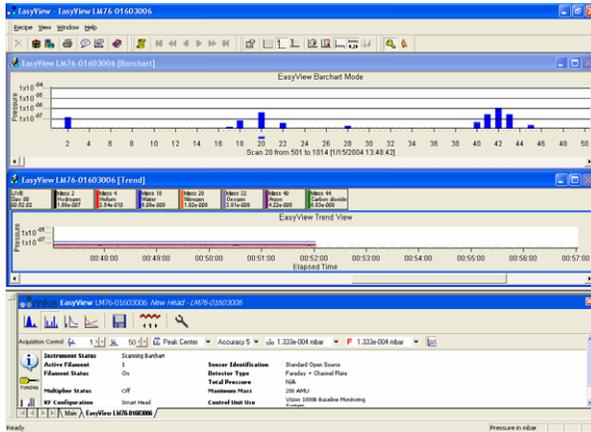

Tile Recipe Vertically

Selecting tile recipe vertically brings the views for the currently active measurements to the front and tiles them vertically.

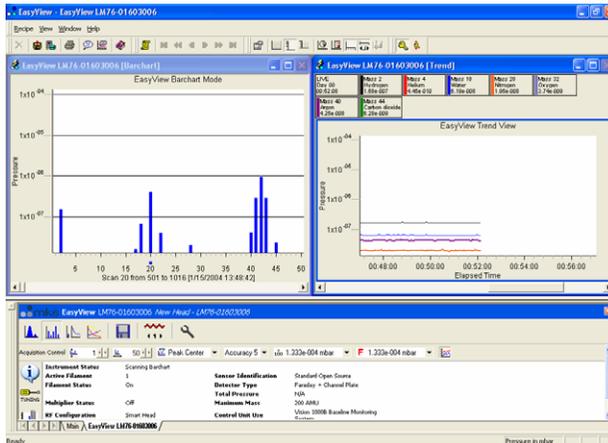

Selecting Individual View

The bottom section of the pull down list contains a list of the currently open views.

The currently active view is highlighted by tick next to it.

To make a different view active, click on that view

9. EasyView Diagnostics

To run the diagnostics mode click on the diagnostics button
The diagnostics mode is only available when no scan mode has been selected.




Running diagnostics produces the following report:

Instrument Diagnostic Information						
15/10/2003 11:41:40						
Serial Number: LM76-00399001						
User Defined Name: New Head - LM76-00399001						
Process Eye Version: 4.00000420						
Instrument Info						
Sensor Identification	Standard Open Source					
Detector Type	Faraday + Channel Plate					
External Hardware	None					
RF Configuration	Smart Head					
Maximum Mass	200 AMU					
Control Unit Use	Standard RGA					
Embedded Software Version	V3.70					
Active Filament	1					
Diagnostics Results						
Diagnostic	Min	Max	Reading	Passed		
-450V Supply	-533.80	-355.90	-409.99	✓		
-15V Supply	-16.50	-13.50	-15.01	✓		
-130V Supply	-143.40	-116.20	-131.80	✓		
+5V Supply	4.75	5.35	4.96	✓		
+15V Supply	13.50	17.50	14.90	✓		
External Pressure	0.00	0.00	0.00	✓		
Source Settings						
	Emission (mA)	Electron Energy (eV)	Ion Energy (eV)	Extract (V)	Faraday A/Inbar	
Standard Electron Energy	1.00	70.00	5.50	-112.00	1.50e-004	
Low Electron Energy	1.00	40.00	5.50	-112.00	1.50e-004	
User defined 1	1.00	70.00	5.50	-112.00	1.50e-004	
User defined 2	1.00	70.00	5.50	-112.00	1.50e-004	
User defined 3	1.00	70.00	5.50	-112.00	1.50e-004	
EasyView	1.00	70.00	5.50	-112.00	1.50e-004	
Mass Alignment/Resolution Settings						
	Low Mass Alignment	Low Mass Resolution	High Mass Alignment	High Mass Resolution		
Standard Electron Energy	32767	32767	32767	32767		
Low Electron Energy	32767	32767	32767	32767		
User defined 1	32767	32767	32767	32767		
User defined 2	32767	32767	32767	32767		
User defined 3	32767	32767	32767	32767		
EasyView	32767	32767	32767	32767		

The report contains information about the control unit configuration, ion-source settings and mass alignment/resolution settings as well as power supply levels. This information can help in fault diagnosis if your RGA is not performing as expected.

In order to test the multiplier power supplies, the filament needs to be switched on. If you run diagnostics without switching the filament on the following dialog appears.

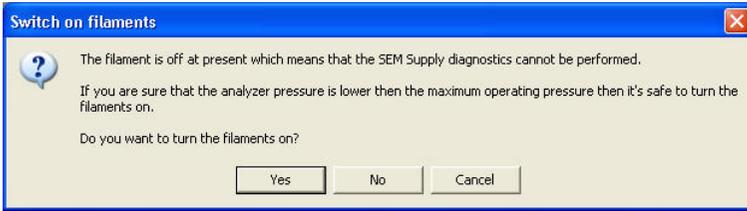

Click <Cancel> to exit without running diagnostics

Click <No> to run diagnostics with the filament off, in which case the multiplier voltage will not be tested.


Click <Yes> to switch the filament on and then run diagnostics.

10. EasyView Tuning

Periodic checks on the tuning accuracy of the instrument should be performed as part of your systems maintenance. Depending on the accuracy of the results you wish to obtain, this period may be between 1 or 12 month intervals.



Note: The tuning parameters can only be changed whilst in the analog mode. In all other modes the tuning information can only be viewed.

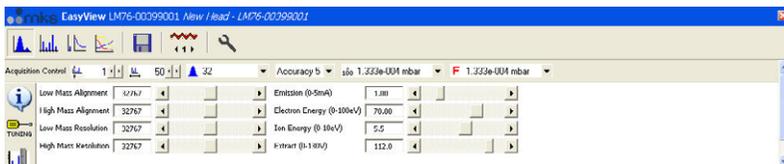

There are two main areas concerning the tuning of the instrument, the mass alignment and resolution controls and the ion-source parameters.

Mass alignment

The mass alignment controls adjust the position of the peaks in relation to the mass scale.

There are two controls for the alignment:

The low mass control affects all masses by the same amount. Moving to slider to the right causes the peaks to move to the right and causes the alignment value to increase. Moving the slider to the left, causes the peaks to move to the left and the alignment number to decrease.

The high mass control affects high masses more than the low masses, for example, moving mass 10 by 0.1 amu using the high mass control, causes mass 100 to move by 1 amu.

Mass resolution

The resolution controls adjust the peak width.


As you increase the resolution the peak will get narrower, but will also become smaller. Reducing the resolution will make the peaks bigger but can lead to interference between adjacent peaks.

The factory settings for standard instruments have the resolution set so the peak width at 5% (10% for HPQ2 units) of the max peak height is 1 amu.

See the following diagram.

This peak is aligned at mass X

The resolution is set to 1 amu at 5% of max peak height.

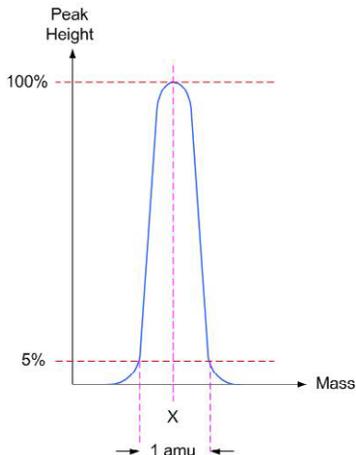


All instruments have the mass alignment and resolution set before leaving the factory. A copy of the tuning numbers (QDAC values) will always be included with the documentation. Only small changes to the alignment and resolution should ever be necessary to account for small changes in temperature between the factory and place of operation. If major changes are made to the settings so that peaks are no longer visible, please re-enter the factory defaults to try and recover the peaks.

11. EasyView Calibration

Mass spectrometers measure an ion-current and not an actual pressure. In order to convert this ion-current to a pressure, we need to calculate a calibration factor.

In calibration mode, the user supplies the actual partial pressure of a given gas and EasyView calculates the calibration factor required to convert the measured ion-current to this pressure.

To view the calibration page click on the calibration button

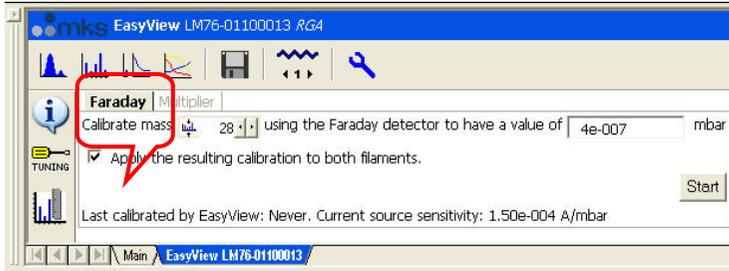


Click on the <Faraday> tab to enter the Faraday calibration setup page.

Choose a mass that you know the actual partial pressure of to calibrate on, and enter the partial pressure of that mass.
To start the calibration click on <Start>


The date of the last calibration using these settings is displayed at the bottom of the window.

Calibration can only be performed if the instrument is not in one of the scanning modes.


The calibration routine will then measure the zero, the peak maximum value and the peak centre value of the mass chosen and then perform a stability check of 10 scans.

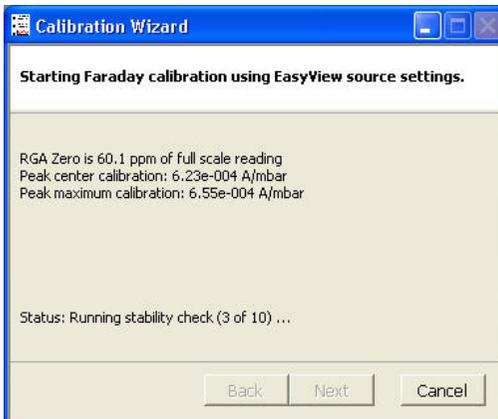


Click <Next> to view the results of the calibration

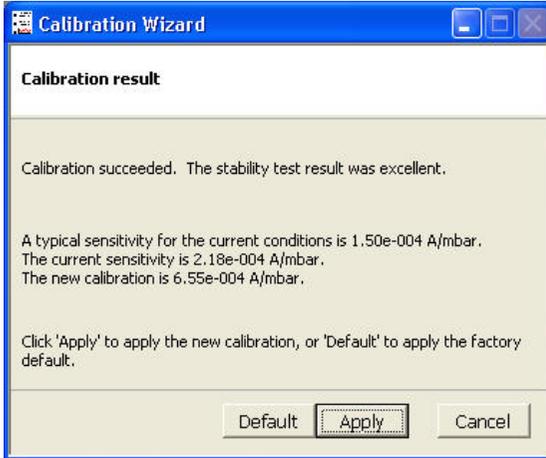

To exit without writing the new calibration information, click <Cancel>. To apply the factory default calibration click <Default>. Otherwise choose <Apply> to write the calibration information.

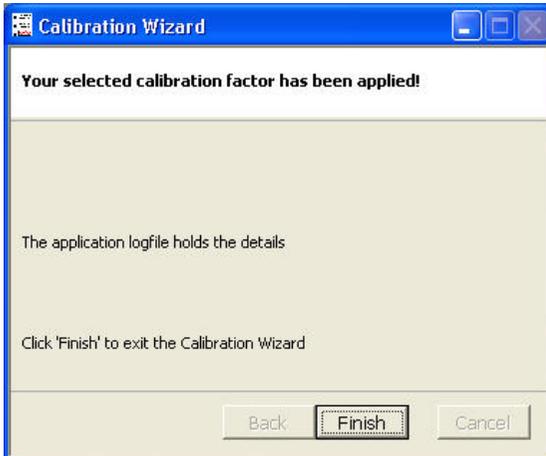

The calibration is now complete click on <Finish> to exit.

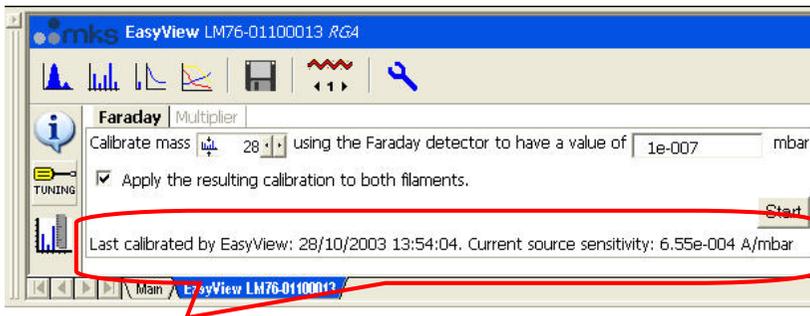

The date and sensitivity value of the calibration just performed is displayed at the bottom of the EasyView window.

```


<MESSAGE Time="2003-10-28 13:52:58" Value="Calibrating RGA ( Source: EasyView, Faraday detector" />
<MESSAGE Time="2003-10-28 13:52:58" Value="Parameters: mass=28;pkht=1.00e-007" />
<MESSAGE Time="2003-10-28 13:52:58" Value="Existing sensitivity: 2.18e-004 A/mbar" />
<MESSAGE Time="2003-10-28 13:53:05" Value="Filaments have been on for 2.8 minutes" />
<MESSAGE Time="2003-10-28 13:53:08" Value="Gain: 100, fsp: 9.16e-005 mbar, RGA Zero is 60.1 ppm of full scale reading" />
<MESSAGE Time="2003-10-28 13:53:10" Value="Reading: 2.91e-007 mbar" />
<MESSAGE Time="2003-10-28 13:53:10" Value="Peak center calibration: 6.23e-004 A/mbar" />
<MESSAGE Time="2003-10-28 13:53:29" Value="Peak maximum calibration: 6.55e-004 A/mbar" />
<MESSAGE Time="2003-10-28 13:53:48" Value="Calibration succeeded. The stability test result was excellent." />
<MESSAGE Time="2003-10-28 13:53:57" Value="Calibration applied: EasyView: New factor = 6.55e-004 A/mbar" />

```

The details of the calibration shown above, are also be logged to the log file, located in `..\EasyView\Logs\Process.xml`.


Multiplier Calibration

To calibrate the multiplier click the <Multiplier> tab on the calibration page.

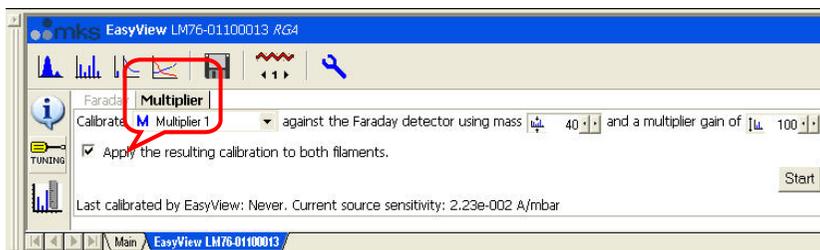


Select the multiplier setting to be calibrated from the drop down list, the mass to calibrate on and the multiplier gain required. The multiplier will be calibrated at the mass chosen, so that the pressure reported in faraday and multiplier at this mass will be the same.


Click <Start> to begin the calibration.

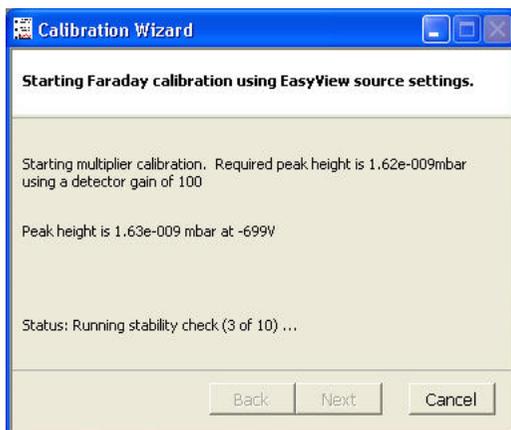


The calibration will do a zero, measure the peak height in faraday, adjust the multiplier voltage so the peak height in multiplier with the gain specified matches and then check the stability.

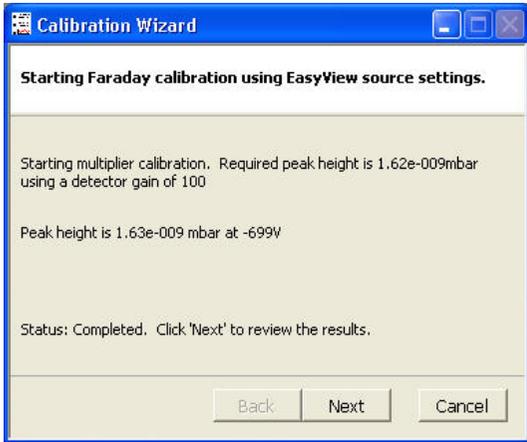

When the calibration finishes, click <Next>.

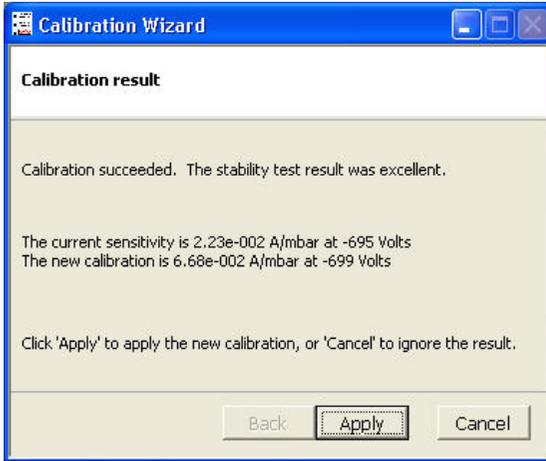

Details of the previous and current calibration will be displayed, click <Apply> to apply the current calibration.

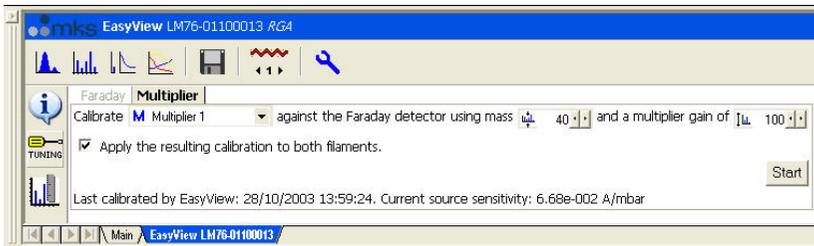


The date and sensitivity of the current calibration will be displayed at the bottom of the EasyView window

```
<MESSAGE Time="2003-10-28 13:52:58" Value="Calibrating RGA (Source: EasyView, Faraday detector)" />
<MESSAGE Time="2003-10-28 13:52:58" Value="Parameters: mass=28;pkht=1.00e-007" />
<MESSAGE Time="2003-10-28 13:52:58" Value="Existing sensitivity: 2.18e-004 A/mbar" />
<MESSAGE Time="2003-10-28 13:53:05" Value="Filaments have been on for 2.8 minutes" />
<MESSAGE Time="2003-10-28 13:53:08" Value="Gain: 100, fsp: 9.16e-005 mbar, RGA Zero is 60.1 ppm of full scale reading" />
<MESSAGE Time="2003-10-28 13:53:10" Value="Reading: 2.91e-007 mbar" />
<MESSAGE Time="2003-10-28 13:53:10" Value="Peak center calibration: 6.23e-004 A/mbar" />
<MESSAGE Time="2003-10-28 13:53:29" Value="Peak maximum calibration: 6.55e-004 A/mbar" />
<MESSAGE Time="2003-10-28 13:53:48" Value="Calibration succeeded. The stability test result was excellent." />
<MESSAGE Time="2003-10-28 13:53:57" Value="Calibration applied: EasyView: New factor = 6.55e-004 A/mbar" />
<MESSAGE Time="2003-10-28 13:55:51" Value="Calibrating RGA (Source: EasyView, Multiplier 1 detector)" />
<MESSAGE Time="2003-10-28 13:55:51" Value="Parameters: mass=40;pkht=0.00e+000;dgain=100;mmpkht=0" />
<MESSAGE Time="2003-10-28 13:55:51" Value="Existing settings: sensitivity: 2.23e-002 A/mbar, voltage: -695.0 volts" />
<MESSAGE Time="2003-10-28 13:56:04" Value="Filaments have been on for 5.8 minutes" />
<MESSAGE Time="2003-10-28 13:57:01" Value="Peak height of mass 40.03amu using Faraday detector: 1.62e-009mbar" />
<MESSAGE Time="2003-10-28 13:57:01" Value="Starting multiplier calibration. Required peak height is 1.62e-009mbar using a detector gain of 100" />
<MESSAGE Time="2003-10-28 13:58:16" Value="The current sensitivity is 2.23e-002 A/mbar at -695 Volts The new calibration is 6.68e-002 A/mbar at -699 Volts" />
<MESSAGE Time="2003-10-28 13:58:32" Value="Calibration applied: EasyView: New factor = 6.68e-002 A/mbar" />
```

The details of the calibration shown above, are also be logged to the log file, located in \.\EasyView\Logs\Process.xml.

12. Preferences

The Preferences menu can be accessed from the <recipe> pull-down menu.

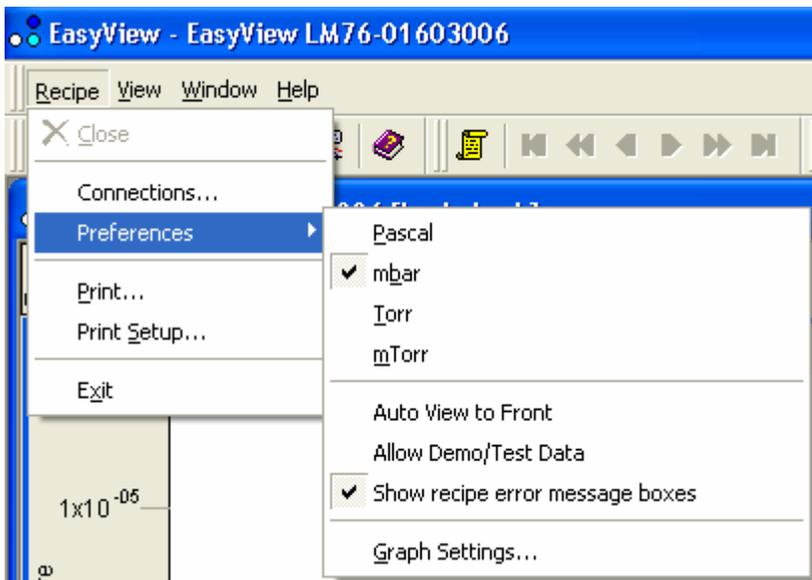

Units of Pressure

EasyView supports the following pressure units, Pascal, mBar, Torr and mTorr.

The internal units used by EasyView are Pascal and all internal pressure values are stored in these units. This enables EasyView and Recall to use different units to display pressures, depending on the preference set in the particular program. Pascal is simply converted to the required unit, before displaying the information.



A "tick" in the selection indicates the currently used pressure unit. All pressure axis will be labelled with the selected unit.

Auto View to front


When EasyView is running a measurement, the view for that measurement is displayed in the Views Window. When more than one recipe is running however, not all views are displayed.

With the <auto view to front> the view for the currently selected instrument is visible.



To select a recipe click on its tab in the EasyView control window

Graph Settings

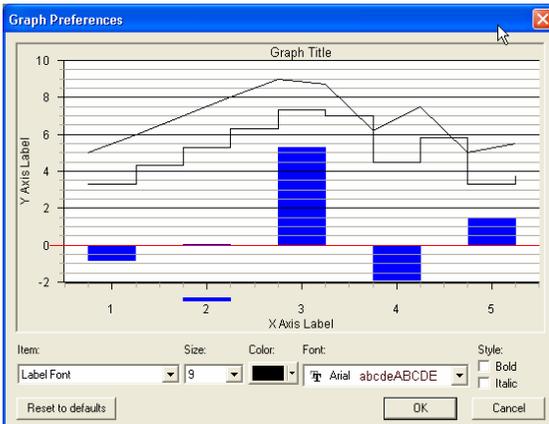

From the <graph settings> option, the global properties for the views can be adjusted.

Common Gases List

The list of common gases which is used throughout the program when editing Trend channels for example, can be tailored to suit your application using Notepad.

This file is located in:

C:/Program Files/Spectra/RGA Applications/Workstation

13. Printing

Both <print setup> and <print> can be accessed from the recipe pull-down menu. Unless a measurement is running, the <print> option will be greyed out. In addition there is a print button on the tool bar.




Printing a View

You may print any active view, if the view is not currently active, click on the view you wish to print, so it becomes the active view. Then either click on the <print> button or select <print> from the recipe pull-down menu.

If you click on the print button, the print will start immediately with the default settings. If you select <print> from the recipe pull-down menu, the standard Windows print dialog appears giving you the option to change the default printer settings.


You may only print swept measurement views and trends, other views within EasyView cannot be printed. If you do require a hard-copy of these other views, use <print screen> from the keyboard.

Below is a printout of a typical Bar Chart spectrum of nitrogen.

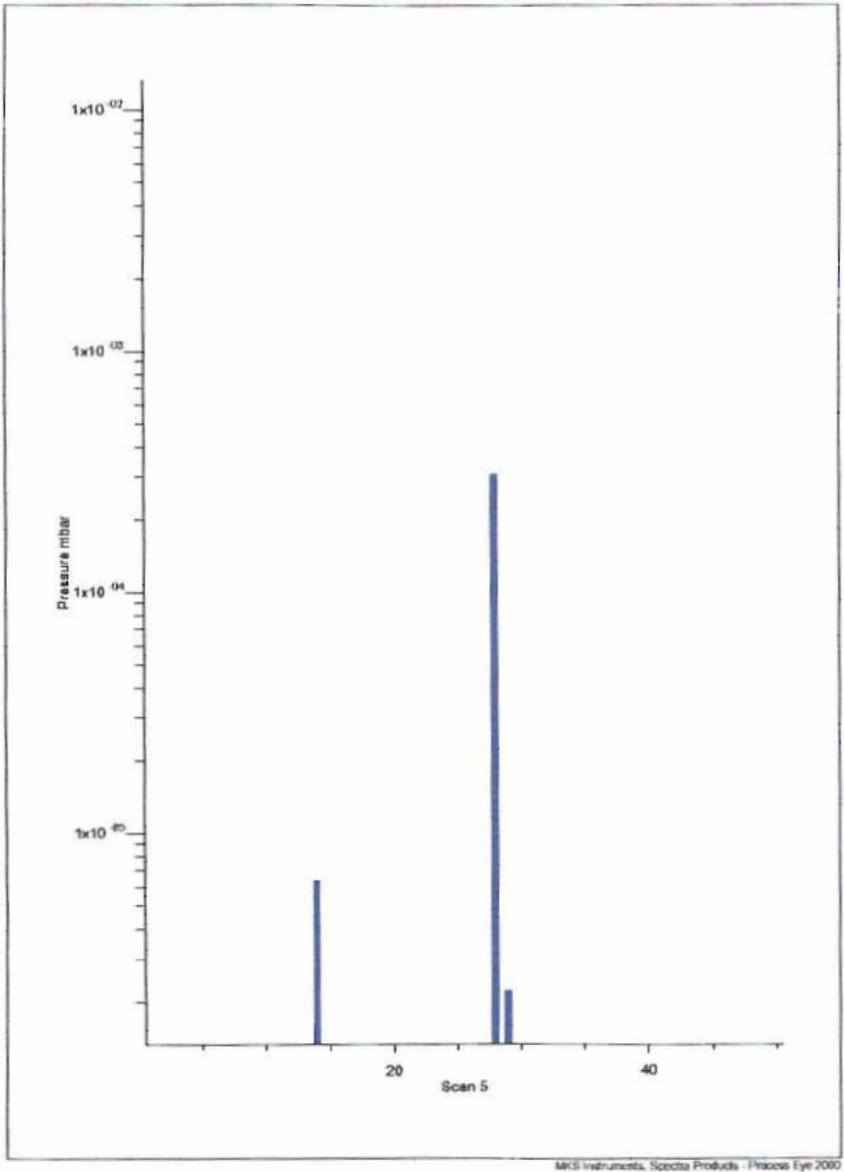

This is a typical trend printout.

The key at the foot of the printout shows the colour of the line, the channel name and the value of the trend line at the start and finish of the trend printout.

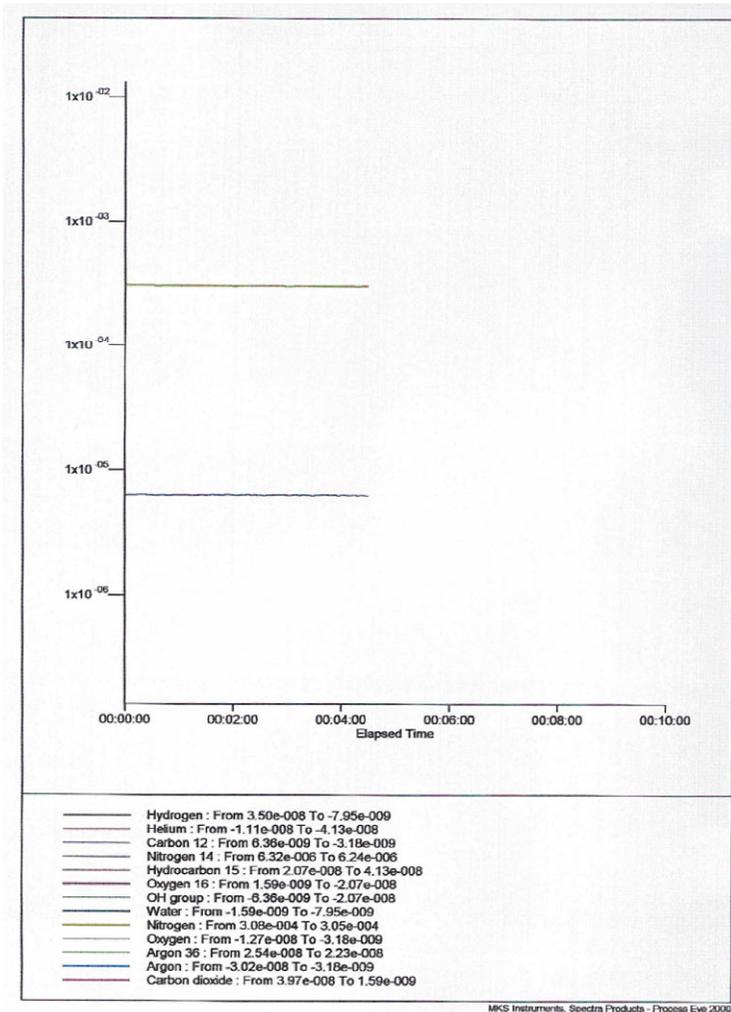

There is no title or foot-note feature in the print facility.

You can however give the printout a title by adding a caption to the view.


14. EasyView Tools

Under the “Tools” menu, there are three programs available as shown. The following section explains each program in turn.



Backup and Recover


The Backup and Recover tool is used to backup the current EasyView installation. The tool will backup the views and registry settings for the current user.

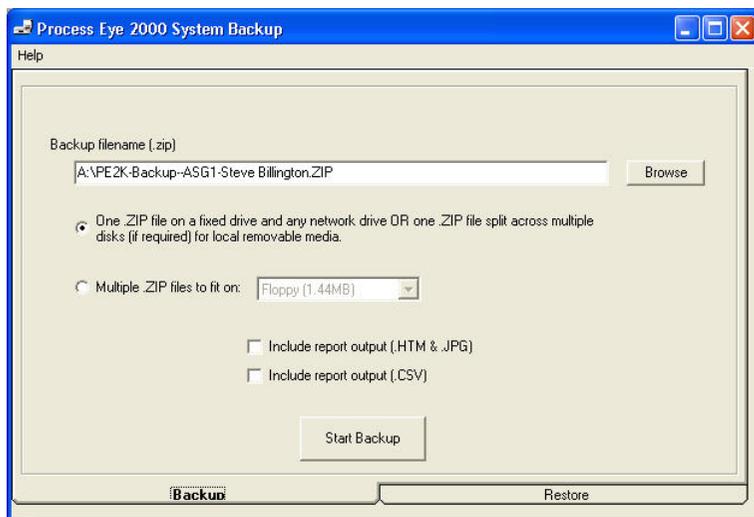

Choose the location and name for the backup file.


To make a spanned backup across a number of disks click the “Multiple Zip files to fit” option and then select the type of media. There is also a custom size option where you may specify the file size of each spanned file.

When ready, click <Start Backup> to begin.

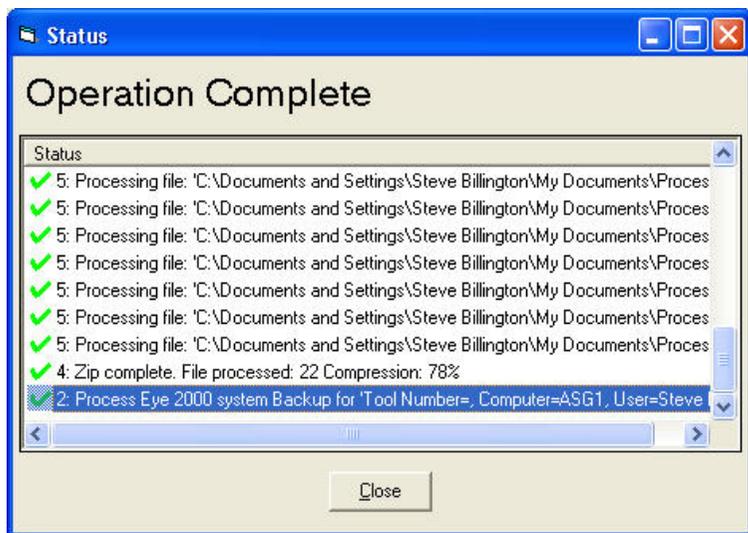


When the operation has completed, click the <Close> button to exit.

To restore from a back-up file, click on the <Restore> tab.

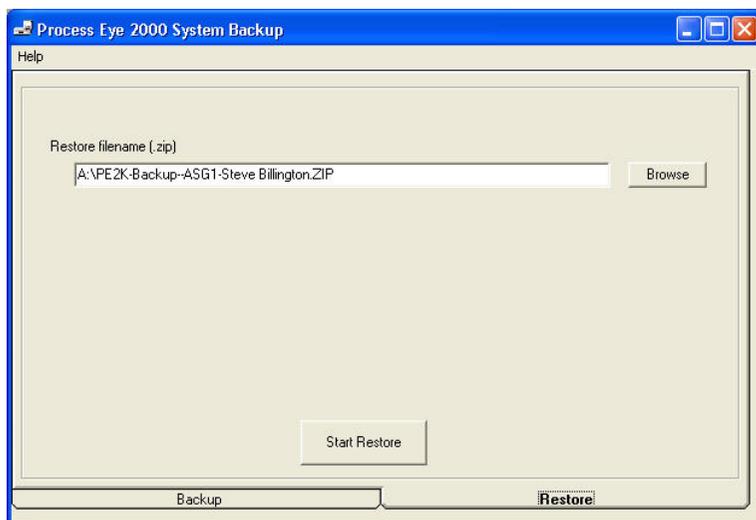

Use the <Browse> button to select the backup and on <Start Restore> to begin.

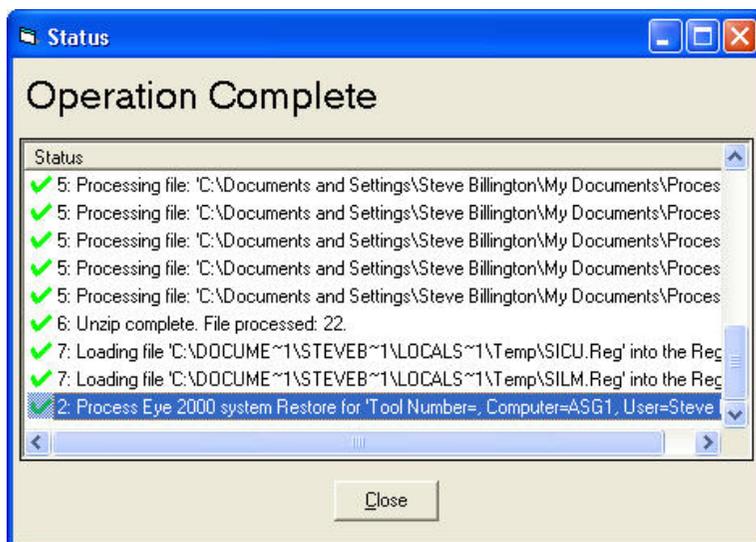


When the restore is complete click the <Close> button.

Computer Lock

EasyView has the ability to lock the computer to prevent unauthorised use.

Note: If you misplace the account information, MKS cannot recover it.

The first time the utility is run a dialog appears detailing the functions of the lock program, click <Yes> to continue.

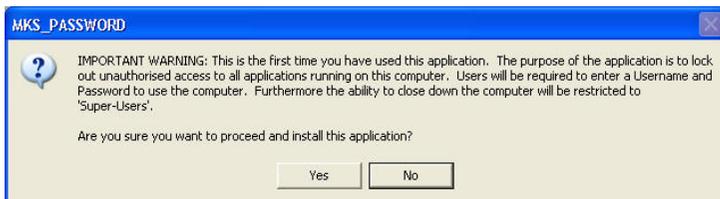

The above dialog displays the initial username of “admin” and a blank password, as the dialog suggests, you should give this “admin” account a new password as soon as possible.

Click <Yes> to continue.



The computer is now locked, to unlock enter “admin” in the UserName text box and then click on <Logon>.




Click on <Configure> to begin adding users, or change the admin password.



In configure you can create, edit and delete users. You can also change the inactivity time before the computer is locked.

Click on <Edit>

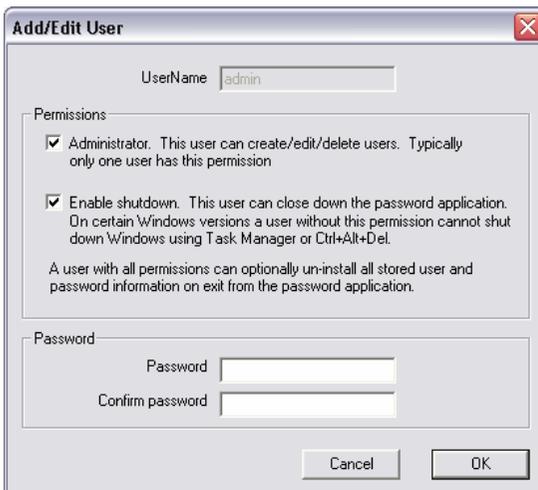

The default "Admin" account should have the options shown selected, if not, select them.

You must always have at least one Admin account with the options shown selected.

Add a password for the admin account.

Click <OK> to save the settings and exit.

To create a new user account, click the <New> button.



Add a "Username", select the permissions and enter a password for the account. Click <OK> to save the account.

Permissions

"Administrator" – can add, edit and delete users.

"Enable Shutdown" – allows the account to close the password application.

Checking both options gives the account the option to not only end the computer lock, but also to either save, or delete the current list of users on exit.

Exiting Computer Lock


Log on using your admin account.



Click on the close button at the top right-hand corner of the dialog.



Click <Yes> to exit the lock, if no user action is taken, the program will revert back to a locked state.

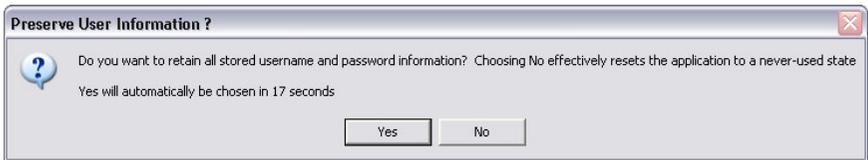

At this dialog, clicking <No> will delete any users, except for the default admin account, you have created.

Clicking <Yes> saves any new accounts you have created.

15. Java Web Applet

It is possible to view and manipulate certain aspects of the EasyView package remotely, via your Internet browser. This provides the facility of network users, without the EasyView software installed, access to the sensor data.


The applet is highly intuitive and includes all the data display routines required for troubleshooting, including leak detection mode.

All that is required is a Java enabled Web browser and the IP address of the unit you wish to view.


Simply enter the IP address of the sensor you wish to view into the "Address" field of your browser and you are taken to the sensor's home page. A Java applet can be downloaded from this page to enable access.

15.1. Accessing the instrument

Start your web browser and type in the IP address of the instrument you wish to view.

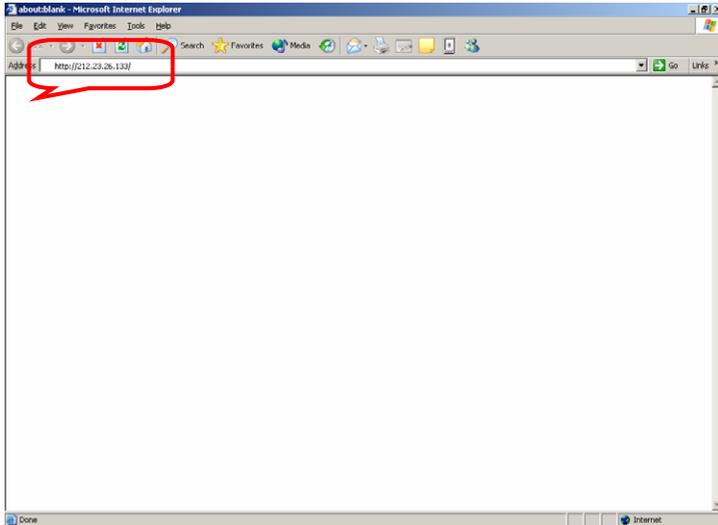


You are directed to the instruments home page.

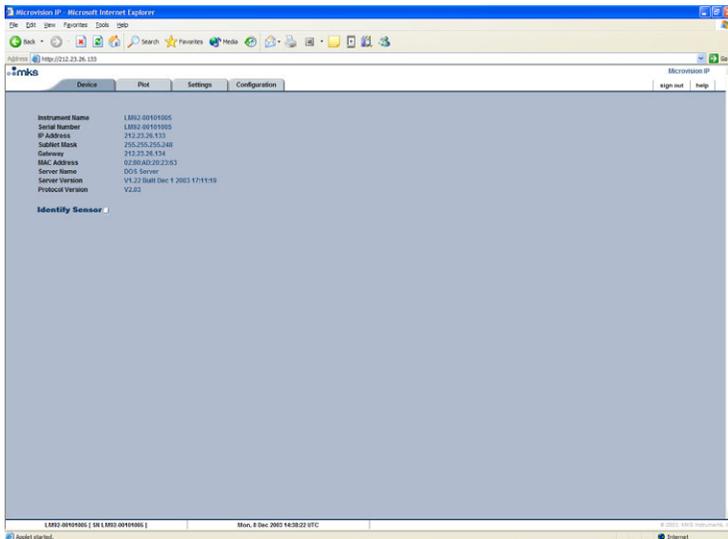


If you chose not to install the Java runtime when you installed the main EasyView program, you will need to do this now before you can proceed.

If a connection to the internet is available, you can download the applet by clicking on the "Plot" tab.

If your PC does not have access to the internet, you will have to install the Java runtime from the EasyView installation disk.

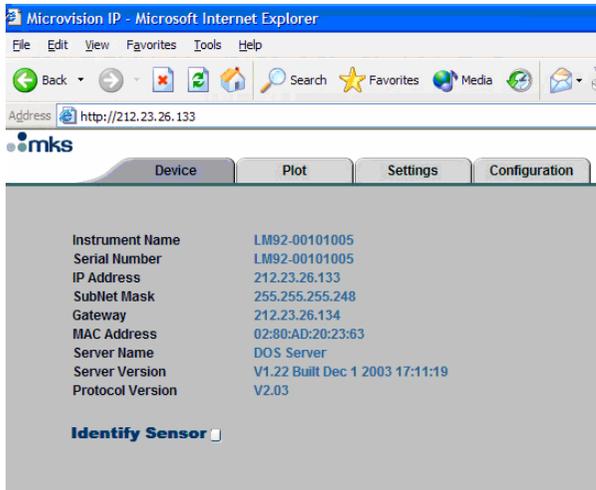


Clicking the "Plot" tab begins the Java download.

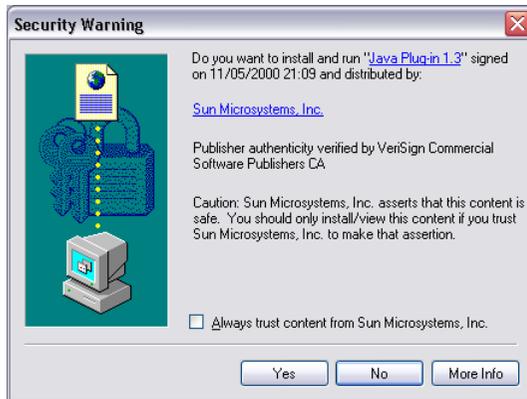

Accept the security warning to begin.


Once the Java runtime has been downloaded and installed, you may need to click the "Refresh" button on your browser.

You now have access to the instrument, to start, click the "Plot" tab once again. The view changes to show the following screen and access to the toolbar is now available.

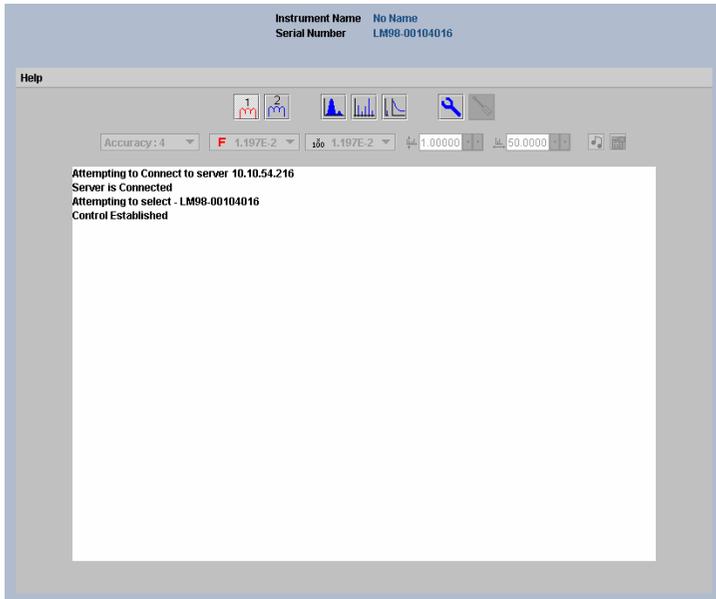


15.2 Page Views and Controls

The Tool Bar

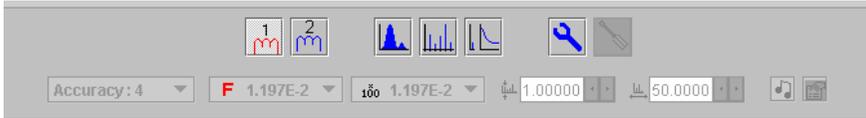


The measurement modes, scan parameters and instrument diagnostics can be accessed from the toolbar. Each function is described in turn through the following section.

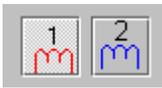

Filament Control

Click the buttons to turn on, or switch between the filaments. The currently active filament is displayed in red.

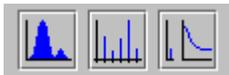

Scan Modes

Analog, Bar Chart and Leak Detect modes.



Diagnostics

Displays a diagnostic report on the instrument.



Tuning


Enter the tuning mode. You must be in Analog Scan mode, this setting is password protected.



Accuracy

Scan accuracy determines the sampling amount carried out on each reading.

A high accuracy number provides a clean scan but increases scan time. The accuracy ranges available are from 0 to 8.

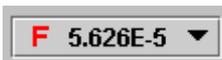


Detector Select

Depending on the instrument configuration, Faraday only, or Faraday and Multiplier detectors are available.




Gain Select

X100 or x20k electronic gain ranges are available.

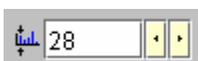

Start Mass

The start mass of the measurement, or when in Leak Detect mode, the probe mass.

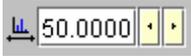

Mass Span

The first mass plus the mass span cannot exceed the maximum mass allowed by the control unit.



Pressure Unit Select

Select from the list of available pressure units.



Tone On/Off

Generates an audio tone in Leak Detect mode. Not available with some sensor packages.

If a more detailed explanation of any of the tool bar buttons is required, these can be found in the corresponding sections of the main EasyView manual.

Page Tabs




Device


Instrument Name	No Name
Serial Number	LM98-00104016
IP Address	10.10.54.216
SubNet Mask	255.255.255.0
Gateway	0.0.0.0
MAC Address	02:80:AD:20:2F:DE
Server Name	Var Not Recognised
Server Version	D1.0 Built Apr 8 2004 13:28:47
Protocol Version	D1.00

Instrument information such as friendly name, IP address and internal software versions are displayed.

Plot

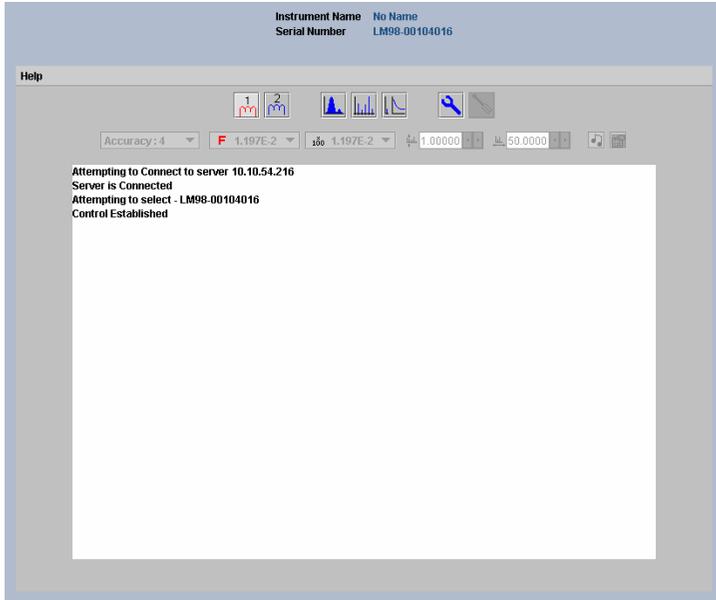

This is the measurement page where most of your time will be spent. From here you can access any of the measurement modes, filament control and diagnostics.

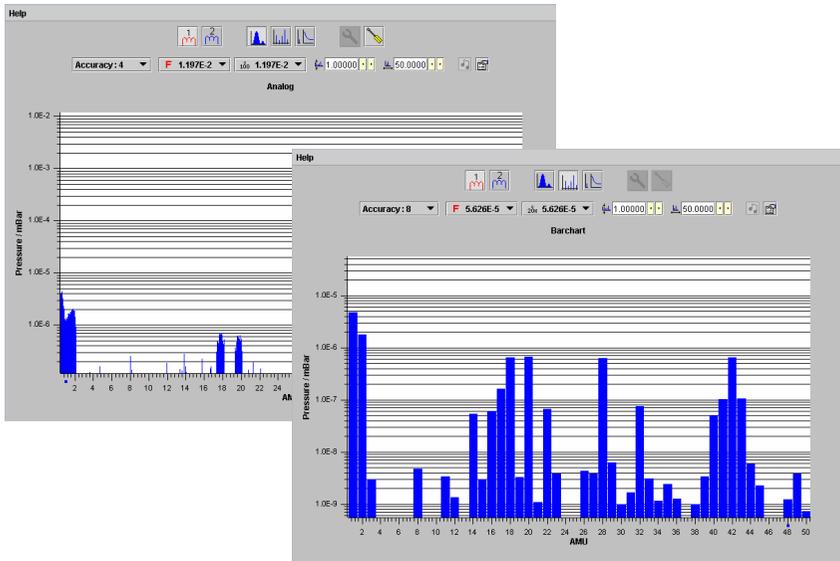

Settings

The Settings page is password protected with a default password of “MKS”. It is recommended that this password be changed to prevent unauthorised changes to the sensor configuration files. See “Security” for details.

**This section is Password Protected.
Please Enter password to continue.**

Enter Password

Sensor **Network** **Security**

Friendly Name

Filament Type

RF Trip

Emission Trip

External Trip

Filament Trip

Filament Control

Done

Under the “Sensor” tab, the following options are available:

Friendly Name

Giving the unit a friendly name is especially useful if you have more than one sensor package installed.

Filament Type

Tungsten or Thoria filament types are available depending on your sensor package option. The filament current drawn by the types of filaments is different, so you should ensure that the correct setting is chosen to match your sensor.

RF Trip – recommended Enabled

If the RF stage of the electronics package runs out of tune, an on-screen alarm is generated.

Emission Trip – recommended Enabled

If a problem is detected with filament emission, the filament is turned off.

External Trip

The filaments can be controlled by an external signal applied via the "X-trip" socket. See the e-Vision hardware manual for signal requirements.

If External Trip is selected, the "Filament Trip" and "Filament Control" options become available.

Filament Trip

Turns the filament off when the X-trip condition is met. The filament must be manually switched on once the trip condition is cleared.

Filament Control

The filament will be turned off for the duration of the X-trip condition, once the condition is cleared, the filament will switch back on automatically.

Configure Sensor

Once the options have been configured, click on this button to write the configuration settings to the sensor.

When finished with this dialog, click the <Done> button to exit the dialog.

Network

The screenshot shows a network configuration window with three tabs: "Sensor", "Network", and "Security". The "Network" tab is active. It contains three radio button options: "Specify IP Address" (selected), "Obtain Address Automatically", and "Set IP Address". Under "Specify IP Address", there are three text input fields: "IP Address" (10.10.54.216), "Sub Netmask" (255.255.255.0), and "Gateway" (0.0.0.0). Under "Obtain Address Automatically", there are three radio buttons: "DHCP", "BOOTP" (selected), and an unlabeled one. A "Done" button is at the bottom right.

Under the "Network" tab, the following options are available:

Specify IP Address

You may change the sensors IP address, Subnet mask and Gateway if required. For help on understanding and assigning IP addresses, please see the on-line help included on the EasyView installation disk.

Obtain Address Automatically

DHCP - Obtain an IP address from your DHCP server.

BOOTP -


Set IP Address

Once the options have been configured, click on this button to write the configuration settings to the sensor.

When finished with this dialog, click the <Done> button to exit the dialog.

Security

The default "Settings" password can be changed here. Enter and confirm your new password.

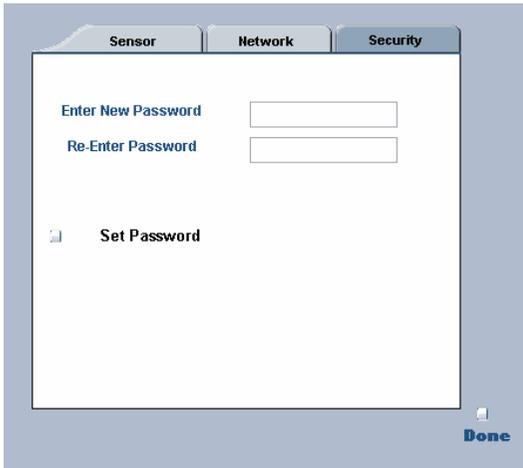

The screenshot shows a software interface with three tabs at the top: "Sensor", "Network", and "Security". The "Security" tab is selected. Below the tabs is a white rectangular area containing two text input fields. The first field is labeled "Enter New Password" and the second is labeled "Re-Enter Password". Below these fields is a button labeled "Set Password" with a small icon to its left. In the bottom right corner of the dialog box, there is a "Done" button with a small icon to its left.

Set Password

Once the password has been changed, click on this button to write the new password to the sensor.

When finished with this dialog, click the <Done> button to exit the dialog.

Configuration

The "Configuration" tab provides a detailed report on the sensor's configuration settings.

The screenshot shows the MKS Configuration interface with the following sections:

- Sensor Information**
 - Serial Number: LM98-00104016
 - Instrument Name: No Name
- General Settings**
 - Mass Range: 100
 - Mass Resolution: Max Res is 1/32
 - Mains Frequency: 50 / 60Hz
 - Detector Type: Faraday Only
 - Filament Type: 0
 - Active Filament: 1
 - X Trip Mode: Trip
 - RF Trip Mode: None Set
- Electronics Settings**
 - Full Scale ADC input: Not Available
 - Full Scale Input Current: 0.000000
 - No. of Electronic Ranges: 2
 - Electronic Gain Ranges: *100 *20k
 - RF Identifier: 1.077881
 - Analog Identifier: -1
- Analyzer Settings**
 - No. of Source Configurations: 1
 - No. of Detector Settings: 0
- Current Settings**

	EE(eV)	IE(eV)	Em(A)	Ex(V)
Standard	-8	8	1	-125
	LMA	HMA	LMR	HMR
Standard	35918	32768	32768	32768

Diagnostics

Serial Number :	Min	Max	Value	Result
-450 Volts	NA	NA	NA	Not Tested
-15 Volts	-16.5	-13.5	-15.195	Passed
-130 Volts	-160.0	-140.0	-149.29	Passed
+5 Volts	4.75	5.35	4.9750	Passed
+15 Volts	13.5	17.5	15.088	Passed
Electron Energy	-68.0	-30.0	-57.739	Passed
Extractor	NA	NA	NA	Not Tested
SEM Supply	NA	NA	NA	Not Tested


Diagnostics can only be performed whilst in the Plot mode without a measurement running.

The diagnostics checks each power supply within the electronics package and displays the result. Some power supplies are not tested, as shown here, depending on the sensor configuration.

16. Communications Troubleshooting

The following flowchart should help you through any connection problems you may encounter.


Note that this assumes you are connecting directly from the host PC to the instrument. If the instrument is to become part of a network, then assistance should also be sought from your IT staff.

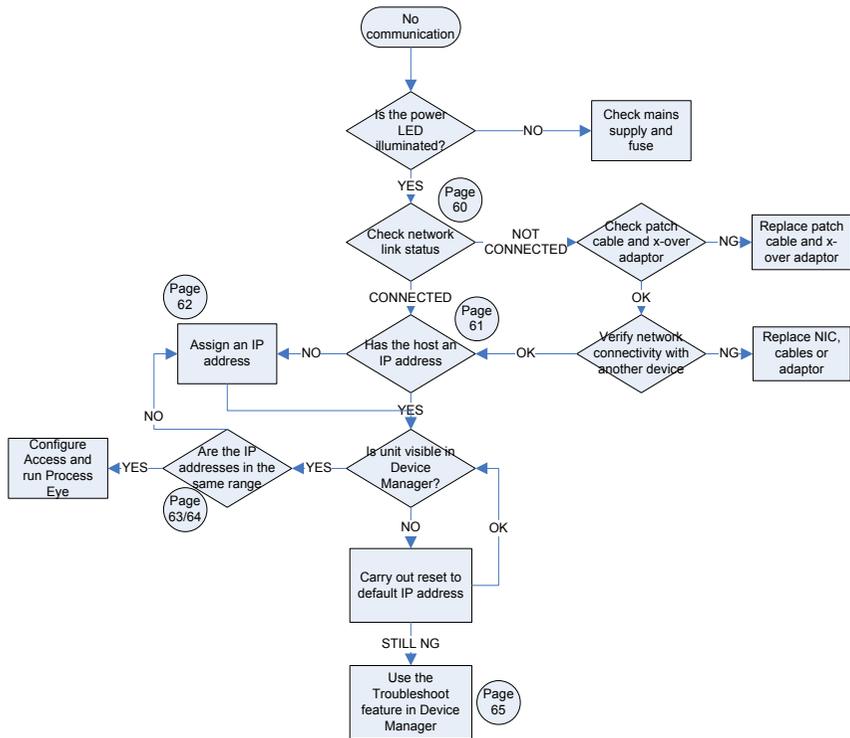

Enabling the Network Connection Notification

To aid in troubleshooting, the network status notification should be enabled. The following paragraph explains the procedure.

Open Windows Control Panel and select the NETWORK CONNECTIONS icon. Choose the network connection you are using for the link with the instrument, right-click and choose PROPERTIES.

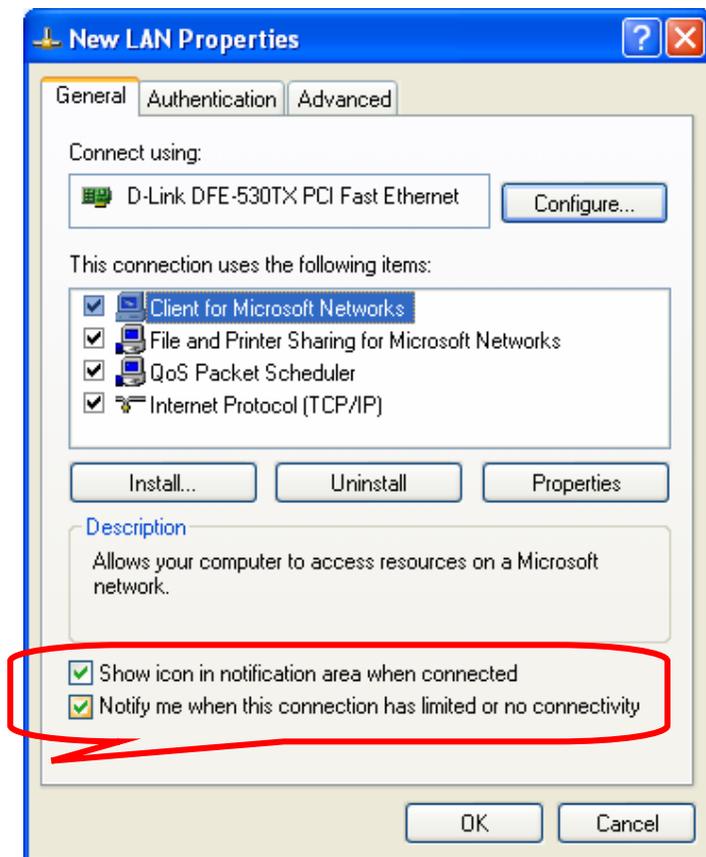

Check the options shown above. Your network connection status is now displayed in the task-bar, near the clock. Mouse-over the icon to check the status of the connection.




Check the IP Address of the Host Computer

Open Windows Control Panel and select the NETWORK CONNECTIONS icon. Choose the network connection you are using for the link with the instrument, right-click and choose PROPERTIES.

Highlight the INTERNET PROTOCOL (TCP/IP) entry from the list and click on the PROPERTIES button.

The following displays a computer configured for automatic IP addressing (DHCP). If you are not part of a DNCP network or intend to directly connect to the instrument, you will need to assign a static IP address.

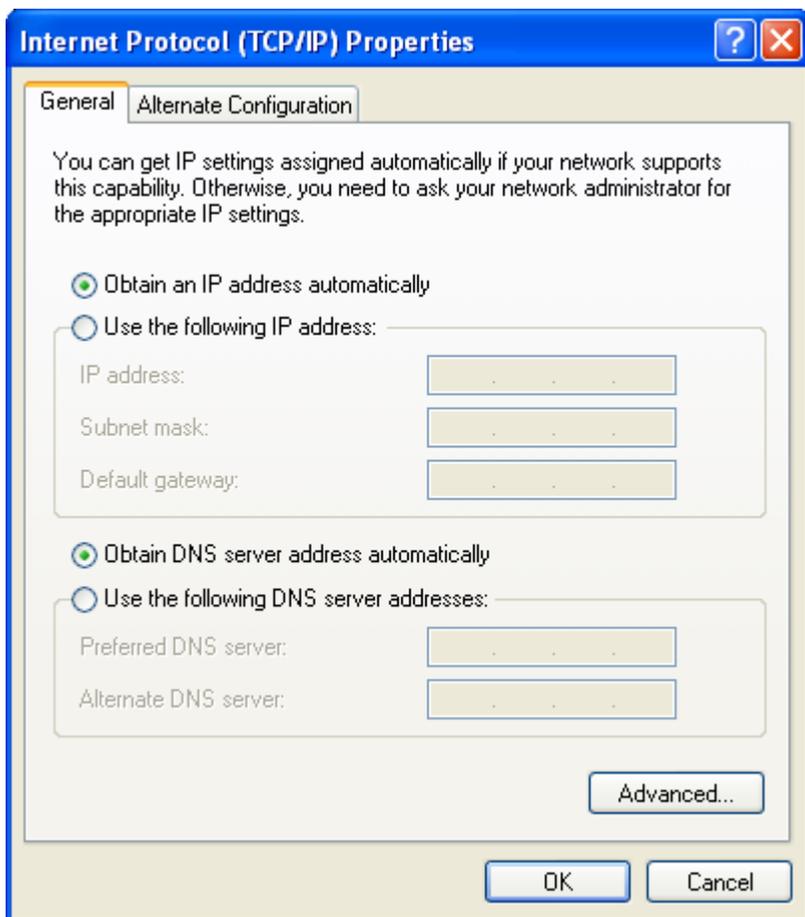


Assign a Static IP Address

Open Windows Control Panel and select the NETWORK CONNECTIONS icon. Choose the network connection you are using for the link with the instrument, right-click and choose PROPERTIES.

Highlight the INTERNET PROTOCOL (TCP/IP) entry from the list and click on the PROPERTIES button.

The default IP address of the instrument is 192.168.0.250, the host PC address needs to be in the same range, but not the same number. As you can see below, this host's IP address ends in 10, but could have been any number up to 255 excluding 250. Use the Subnet shown.

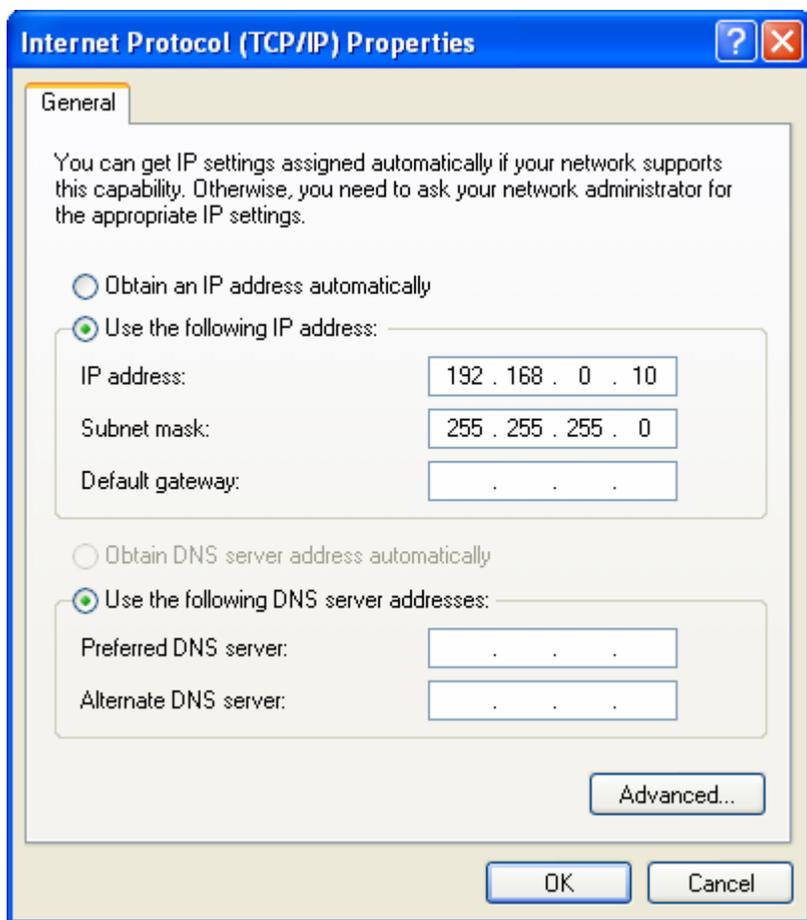


PC and Instrument on different IP ranges

You can use Process Eye's Device Manager to check on the current IP address status of both the host PC and the instrument.

Start DEVICE MANGER, highlight the instrument, click the CONFIGURE button and choose CHANGE THE IP ADDRESS from the list.

You can see in the dialog below, that the IP address of the instrument differs from the IP address of the host computer.

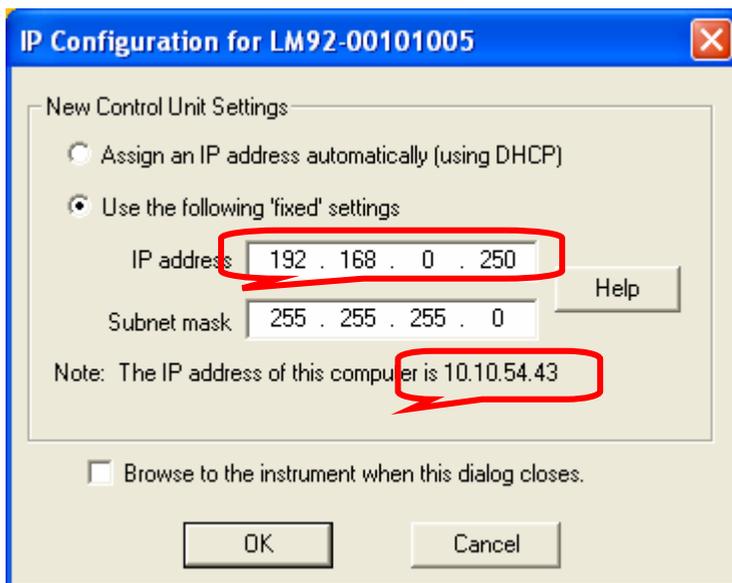

In this example, the host PC will not be able to communicate with the instrument while the two IP addresses are in different ranges.


There are two options. Change the IP address of the host PC, or change the IP address of the instrument.

To change the IP address of the instrument, overwrite the current address with the new one.


Note: If the host PC is part of your network, then changing its IP address is not recommended without seeking advice from your IT staff. Likewise, assigning a new IP address to the instrument must be done under advice or network conflicts could occur.

PC and Instrument on the same IP ranges

You can see in the dialog below, that the IP address of the instrument and the IP address of the host computer are in the same IP range.

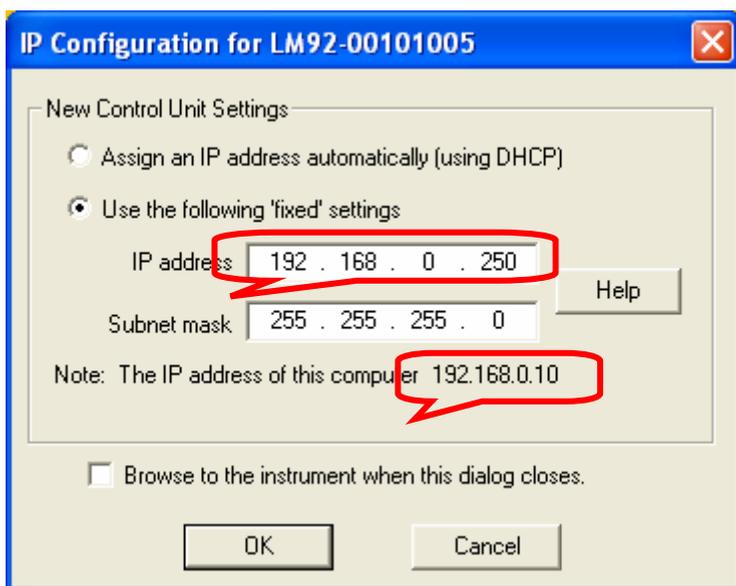

In the above case communication between the host PC and the instrument is possible, though Device Manager will recommend you change the instruments IP address to another, not ending in .250. This is to avoid potential problems when installing more than one instrument.

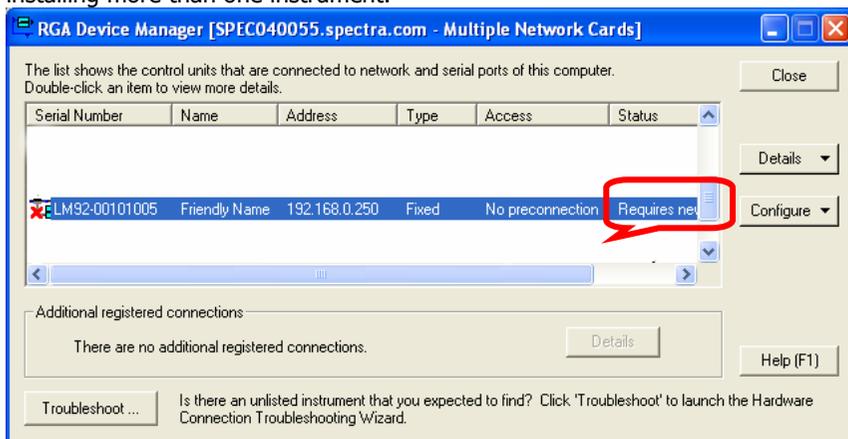

Resetting the Default IP Address

To reset the instrument to its default IP address of 192.168.0.250, carry out the following procedure:

Power on the unit

Press the RESET button and immediately press and hold the MODE button for 20 seconds.

Important:

If the instrument has been configured for DHCP, keep the MODE button pressed for 90 seconds, or a reset will not occur.