




# HPQ2-T

## Additional Setup Instructions


## Installing Process Eye Professional

Carry out the Administrative installation as usual, when you perform the “User Setup” ensure you choose the option to install the “HPQ2-T Getting Started” recipes. This will ensure you have the HPQ2-T Status dialog visible on the default recipes.

If you don't see the option for the HPQ2-T recipes, it is likely you already have Process Eye Professional installed on your machine. In this case, you will need to click the “Have Disk” button as shown.

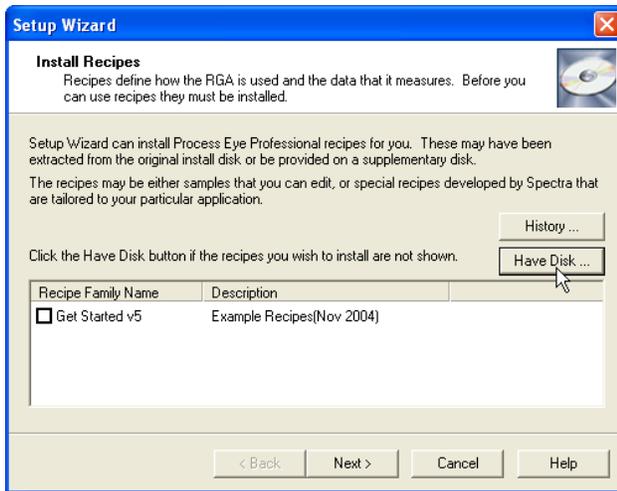


From the pull-down menu choose the “Browse” option

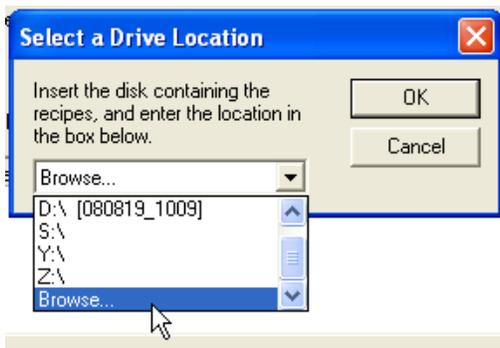


Browse to:

CD drive / Software / LocalMasters / HPQ2T Getting Started Recipes and select the recipe list as shown.

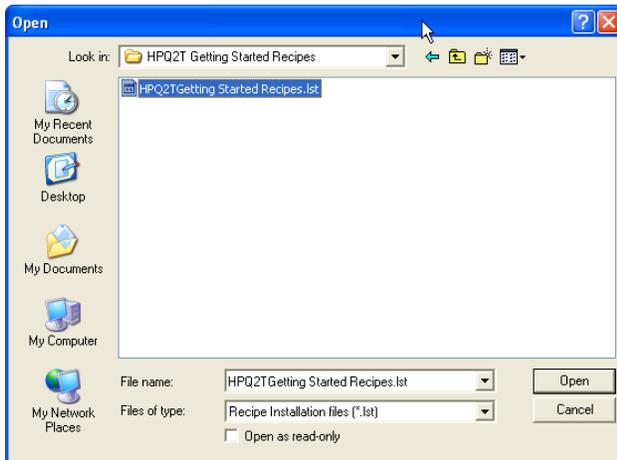

You will now see the HPQ2T recipes are available for selection

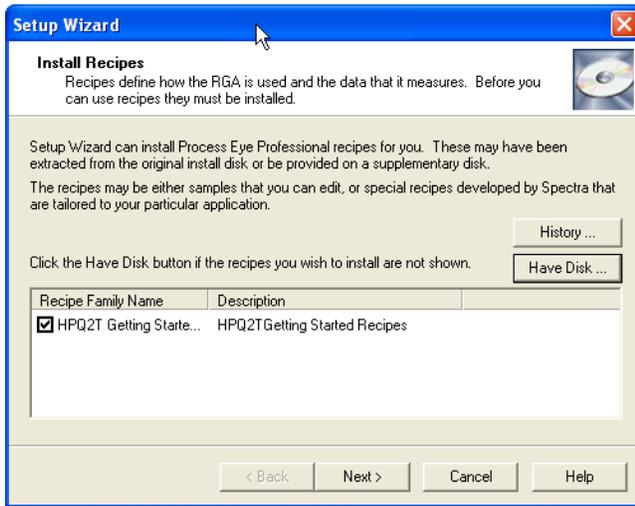


Check the box and click "Next" to continue.

Once Process Eye Professional has been correctly installed there are some settings unique to the HPQ2-T system that will need to be defined:

## *Inlet Factor*

The Inlet Factor can be found on the system information paperwork supplied. Make a note of this value and browse to the following location on the RGA host PC:

If a common location for the Process Eye files option was chosen during the Process Eye installation:

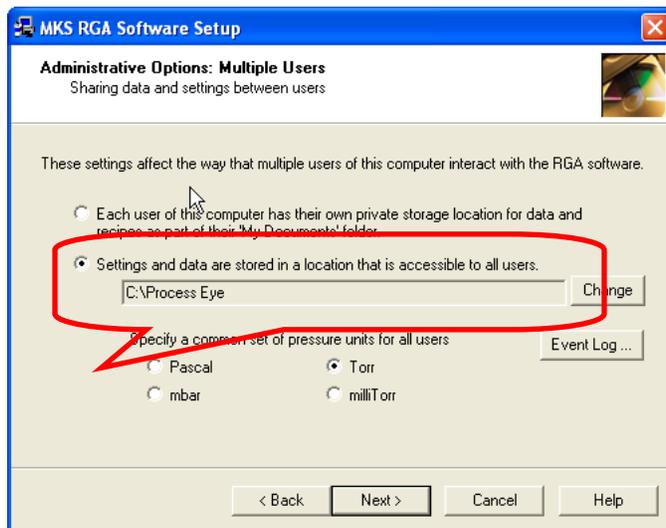

Browse to:

C:/ Process Eye / Views


Open the file named "InletFactors.xml" and look for the following line:

```
<rga name= "Friendly Name" InletFactor= "80" />
```

Change the "InletFactor= "80" to the inlet factor value you noted from the system information paperwork

The "Friendly Name" you intent to call the unit must also be stated here, for ease we suggest "HPQ2T", but this can be any name you wish. However, it must be exactly the same as the friendly name given when setting up Process Eye Professional.

If the option to make files private to each user was chosen during the Process Eye installation:

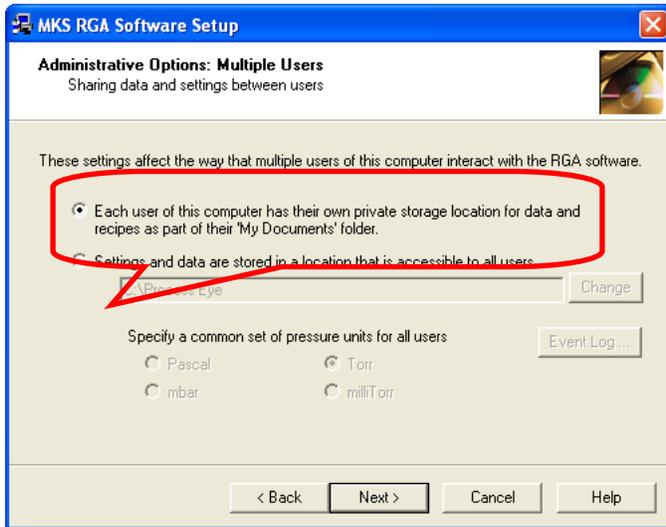

Browse to:

C:/ My Documents / Process Eye / Views

Open the file named "InletFactors.xml" and look for the following line:


```
<rga name= "Friendly Name" InletFactor= "80" />
```

Change the "InletFactor= "80" to the inlet factor value you noted from the system information paperwork

The "Friendly Name" you intent to call the unit must also be stated here, for ease we suggest "HPQ2T", but this can be any name you wish. However, it must be exactly the same as the friendly name given when setting up Process Eye Professional.

## Writing Custom Recipes

If you wish the HPQ2-T Status dialog to be displayed when writing custom recipes, be sure to check the option to use an "Include File" as shown below.

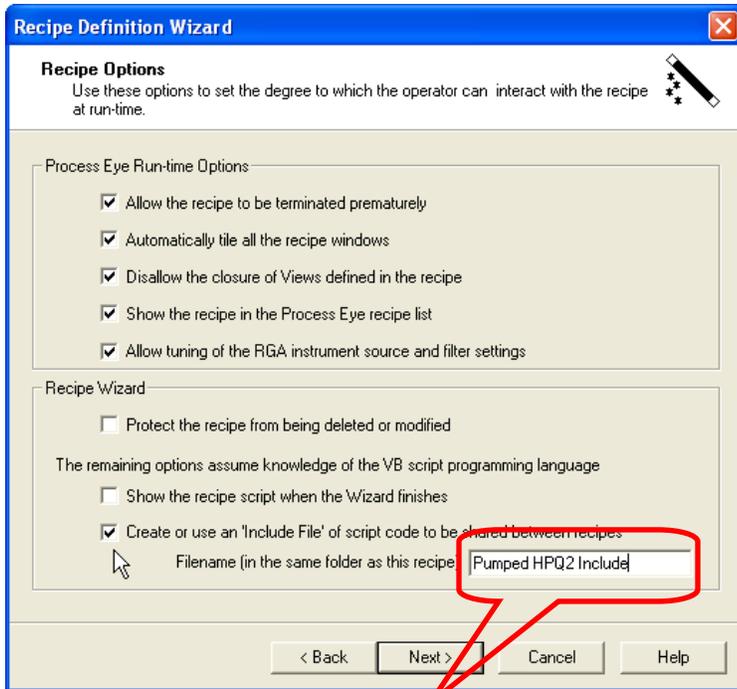

Specify the file name "Pumped HPQ2 Include" then continue on with creating your recipe.