

## ISS-2 Integrated Sampling System

The **ISS-2 INTEGRATED SAMPLING SYSTEM** is a fully integrated 1-cm cuvette holder and tungsten halogen light source. It couples to our spectrometers with optical fiber to create a small-footprint systems for VIS-NIR (~360-1100 nm). The ISS-2 has a 900 hour bulb. Also, in the ISS-2, the fan is not enclosed in the base; it is exposed and should be handled with care.

### Parts Included

- ◆ ISS-2 cuvette holder and light source assembly
- ◆ 12VDC wall transformer for providing power
- ◆ 1-cm square, plastic cuvette for holding samples
- ◆ Screwdriver for adjusting the fit of the cuvette
- ◆ Allen wrench for adjusting the collimating lens

### Operation

#### Attaching the Fibers

1. Attach a SMA-terminated optical fiber to the collimating lens.
2. Attach the other end of this fiber -- the read fiber -- to the spectrometer.

#### Adjusting the Fit of the Cuvette

The ISS-2 is preset for use with a 1-cm cuvette. When properly adjusted, the cuvette should fit snugly into the holder.

1. Locate the two ball plunger screws.
2. Use the screwdriver to loosen the two ball plunger screws until the ends are visible in the holder.
3. Insert your cuvette into the holder.
4. Tighten the screws until the ball contacts the cuvette and starts to compress. Do not over-tighten.

#### Installing Filters

1. Loosen the filter clamping screw with an Allen wrench.
2. Insert the filter into the filter slot. The filter slot can accommodate filters up to 6 mm thick.
3. Clamp the filter in place by gently tightening with an Allen wrench.

#### Turning On the Lamp

1. Plug the wall transformer end into a 110 V outlet. Plug the 12 V output end into the back of your ISS-2.
2. Screw a fiber into the SMA connector.
3. Find the on/off switch on the back of the lamp and turn the lamp on.

#### Replacing the Bulb on the ISS-2

1. Order a replacement bulb.
2. Turn off the lamp and allow it to cool.
3. Remove the fan from the bottom of the lamp by loosening the four screws securing it to the base. Two of the four screws also hold the two front legs in place. Once the screws are loosened, gently pull the fan, along with the legs, away from the lamp.
4. Now that the fan is removed, use an Allen wrench to loosen the set screw that is underneath the fan. This screw holds the bulb in place. You do not need to remove the set screw -- loosening it is sufficient.
5. Locate the two set screws at the back of the lamp, one above each back leg. These two screws keep the two halves of the lamp together. Remove the two screws.

6. Gently separate the two halves of the lamp and pull the bulb out of its housing.
7. Detach the wire and socket from the lamp leads. Remove bulb unit and discard.
8. Plug the new bulb into the socket and slide it forward into the front of the lamp as far as it will go.
9. Tighten the set screw on the bottom of the lamp to hold the bulb in place.
10. Close the two halves of the lamp, being careful not to pinch the wires. Replace the two screws at the back of the lamp and secure the fan to the bottom of the lamp.

## Specifications

Path length:	1 cm
Collimating lens:	BK 7 glass (~360 nm - 2 μm*), 5 mm diameter, f/2
Collimating lens termination:	SMA 905
Filter slot:	accepts filters up to ¼" (6 mm) in thickness
Base material:	aluminum
Spectral range:	360 nm - 2 μm*
Dimensions:	9.0 cm x 5.0 cm x 3.2 cm (LWH) 3.5" x 2.0" x 1.25" (LWH)
Power input:	12 VDC/800 mA, 7-20 VDC/0.5-2 amps
Power output:	6.5 watts
Bulb life:	900 hours
Bulb color temperature:	3100K
Output to bulb:	5 volts/1.3 amps
Output regulation:	0.2% voltage
Time to stabilized output:	~ 30 minutes
Bulb output:	7400 foot-candles (7.4MSCP)

\* Though the product can be used to 2 μm, it can be configured to "see" only to 1100 nm with our S2000 spectrometer.