

FIB-SEM Nanofabrication: Key Strengths for Benchmark Applications

Introduction

Over the last decades, focused ion beam (FIB) technology has evolved from a niche method into an extremely versatile technique for various application areas. FIB's great success has been accelerated in particular by the combination with scanning electron microscopy (SEM) in FIB-SEM microscopes. Today these FIB tools are found in most microscopy and analysis labs worldwide, where the FIB column is included as an add-on component for supporting sample preparation and basic patterning tasks. However, FIB is also an increasingly valuable tool for nanofabrication and rapid prototyping applications, providing direct resistless three-dimensional patterning and thus complementing other top-down lithography instrumentation. Even though patterning with ions has advanced very quickly to a high performance level, the SEM prioritized microscope architecture has limited it so far to more sophisticated applications. Based on unique FIB and precision stage technology as well as a novel nanofabrication platform, the Raith FIB-SEM is designed around the FIB as the primary patterning technique, supported by uncompromised SEM and laser interferometer stage capabilities. Such a FIB-centric state-of-the-art nanofabrication system allows scientists to take advantage of the unique capabilities afforded by focused ions. This paper discusses various capabilities and key strengths, based on exemplary applications from different nanopatterning areas.

Strengths and limitations of EBL

Today's state-of-the-art nanofabrication systems must deliver flexibility, stability over space and time, accuracy, and ease of use. These demands are fulfilled by EBL patterning: precise laser interferometer stage (LIS), best positional stability, enhanced beam current stability, excellent beam size deflection and accuracy are mandatory for these instruments. Moreover, software with industry standard GDSII designs, advanced beam control, navigation, and true automation are essential components of any dedicated nanofabrication system. These attributes have made EBL the most prominent fabrication method in the field of nanofabrication, achieving wafer scale patterning to nanometer accuracy in terms of placement and feature size in a reasonable time. However, EBL unfolds its strengths only in combination with other techniques used in the process chain. For transforming a substrate

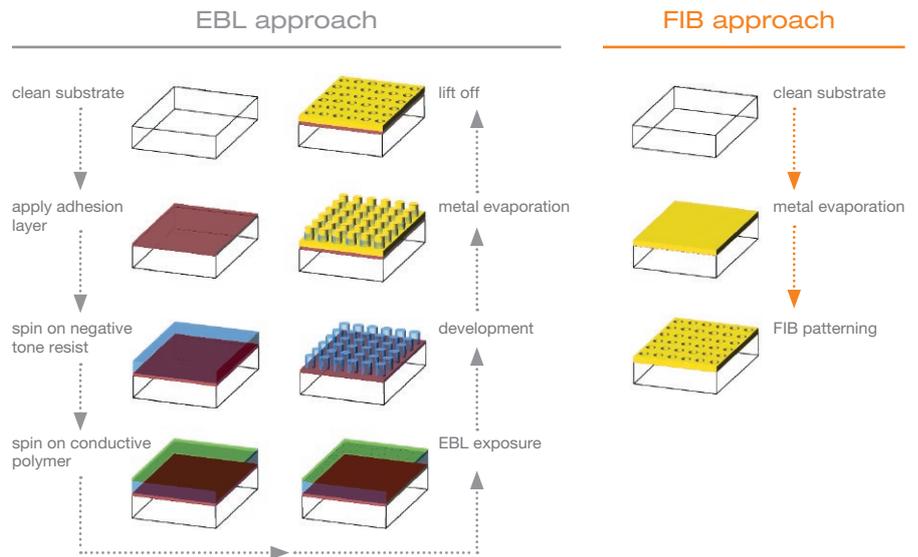


Figure 1: Number of steps for e.g. lift-off process is strongly reduced in case of FIB (right) compared to EBL approach (left).

into a patterned substrate, EBL patterning is just one process step among many others, such as etching and lift-off. Furthermore, EBL is a

superior method for lithography on flat substrates, where patterning of topographic samples can be complicated to achieve.

Lean Nanofabrication: Core process steps addressed by FIB

The limitations of long process chain and patterning of topographic surfaces can easily be overcome by using a focused ion beam to structure substrates. FIB-SEM instruments are commonly used for 3 dimensional patterning – even over topographic samples. As another benefit of these systems, the resistless fabrication method allows direct patterning with far fewer process steps. As a consequence, development of new patterning processes can be achieved very fast thanks to *in-situ* process control.

Conventional FIB-SEM systems, designed for various microscopy and analysis applications in a SEM-centric setup, exhibit some limitations for nanofabrication applications. The FIB column is inclined with respect to the horizontal stage motion and the column and stage technology themselves, and this characteristic, in combination with the overall system and software platform, does not support advanced FIB nanopatterning. Long-term highest-precision automated unattended or large-area processes demand a different class of technology and system architecture. These FIB capabilities plus the *in-situ* SEM process control unlock new pathways and help achieve results more rapidly for an increasing number of applications.

The nanopatterning optimized FIB-SEM system

To advance FIB techniques to the performance level of a lithography tool, Raith's FIB-SEM systems follow a different approach in system architecture. Here, the FIB column is designed as the primary technique. Within a dedicated FIB-centric architecture, it is oriented vertically to the stage, while the SEM column is inclined with respect to the horizontal stage. In addition, the on-board LIS allows highly precise movement of the sample with respect to the ion beam and to the electron beam (see Figure 2).

Operating with Raith's mature nanoFIB Three column, this system concept is the first to offer access to a wide range

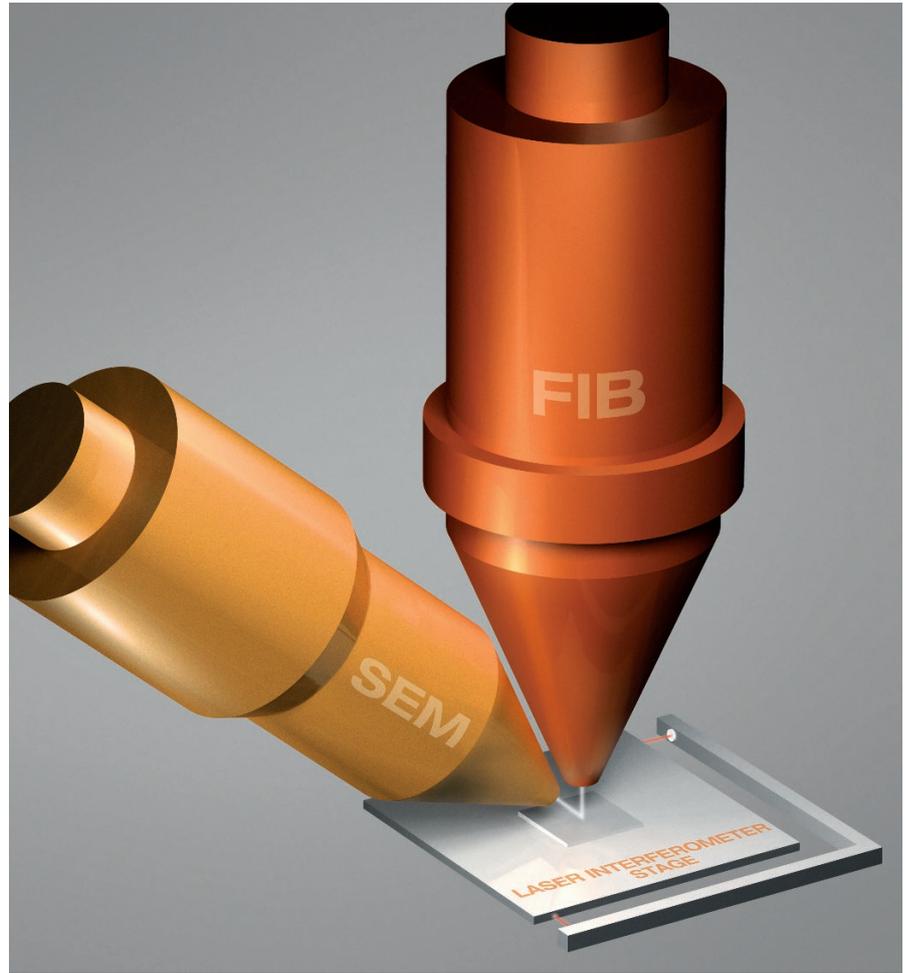


Figure 2: Schematics of Raith's FIB-SEM setup including the highest precision laser interferometer stage

of applications which are beyond the reach of common FIB-SEM instruments:

- Fabrication of complex patterns spreading over multiple deflection fields: this application requires high precision stitching capability of the LIS and, at the same time, extremely low beam distortion within each single write field.
- Furthermore, the ion column is optimized for highest resolution and excellent beam characteristics offering the tightest beam spot of any gallium FIB by far. Only this level of beam performance allows patterning of e.g. dense structures with nanometer gaps in between.

- Wafer scale FIB nanofabrication is even more advanced. In this case, the above mentioned properties are complemented by continuous focus control, so that minimum feature size and optimum placement accuracy are achievable at any position on the wafer.
- Advanced process control of FIB processes can be easily achieved by various *in-situ* techniques (e.g. end-point-detection and SEM imaging).

This approach makes Raith's focused ion beam systems the first truly dedicated FIB-SEM nanofabrication systems. In the following, typical applications illustrate the versatility and high performance of this approach.

FIB hard masking of large area patterns

The first example shows a novel way of employing specific ion beam effects for the purpose of fast patterning. When patterning a substrate using a low ion dose, the Ga ions are implanted into the surface without even milling the top surface layer. Instead of creating topographic structures, the sample surface can be functionalized in a very controlled way by this Ga implantation with high lateral resolution. In a subsequent etching step, this functionalized top layer serves directly as a hard mask, resulting in a very simple and fast process. An example of the process workflow is illustrated in Figure 3.

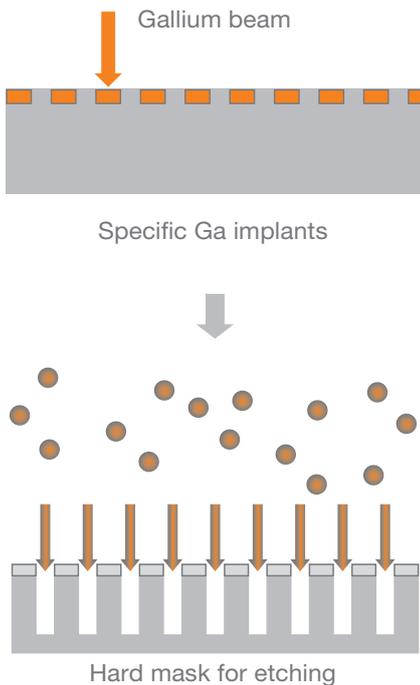


Figure 3: Sketch of the functionalization process – after controlled Ga implantation trenches are etched into the substrate.

After functionalization of a diamond surface with Ga ions and subsequent fluorine plasma etching, a grating with about 400 nm line width and 50% duty cycle was fabricated (see Figure 4). The complete design consists of multiple gratings with varying orientation. The overall structure extends over an area of 3 x 4 mm². The resulting pattern forms a diamond-like structure illustrated in the photograph of Figure 5. The different

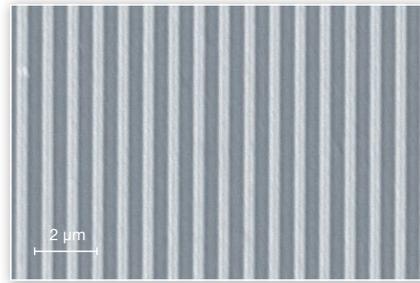


Figure 4: SEM viewgraph of the etched grating [1]

grating orientations generate different colors in the optical viewgraph.

The functionalization step took 4h, which is comparatively fast.



Figure 5: Photograph of the complete pattern after etching. Colors represent the different grating orientations. [1]

Nevertheless, the process requires high stability in terms of beam current, position drift of the ion beam, and laser stage, as well as perfect stitching accuracy. (Any stitching error, however small, would be magnified in the optical viewgraph.) In addition, focus control on the fly is mandatory for achieving perfect structuring conditions over the complete area of the pattern. The example shows how processes which are typically quite complex can be simplified by turning a contamination into a desired functionalization and making use of the outstanding performance of the FIB system, requiring only the substrate and the Ga beam.

High resolution features in a large array

A FIB spot is usually not described by a pure Gaussian distribution, but instead exhibits significant beam tails outside the central spot. The contribution of these low dose beam tails becomes important for gas assisted processes or creates collateral damage close to, or even inside, the desired pattern. Although this dose is very small (10⁻² – 10⁻⁵ of peak intensity), it can still have a significant influence on the result: the superposition of the beam tails at the edge of a pattern or between adjacent and dense patterns can lead to unwanted edge rounding or sample damage. Challenges for achieving small feature sizes and at the same time a large array for final measurements are given e.g. in the field of

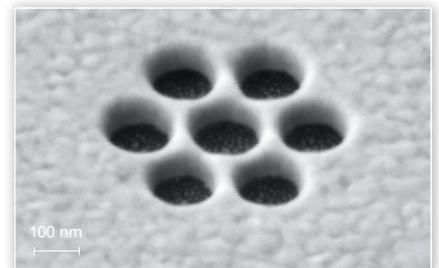
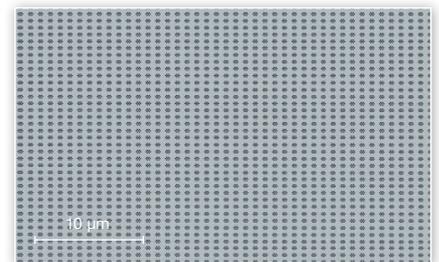


Figure 6: Oligomer structures within a 100-μm-wide array for plasmonics, showing excellent fidelity and reproducibility [2]

plasmonics, where large numbers of identical structures are designed with high pattern density. Figure 6 shows an array of 12,000 oligomers each consisting of 7 circles. Between two adjacent circles, the remaining strip of 80-nm-thick gold has a width of only 30 nm. The ability to create such a strip in gold is also good proof of excellent beam spot characteristics, specifically the record low beam tails, since the beam profile overlaps from several directions at this point. Finally, it is noted that the film consists

of polycrystalline gold, which is even more difficult to pattern than single crystal gold films in terms of high resolution and reproducibility.

Development of an advanced FIB process

Fresnel zone plates (FZP) for X-rays require high accuracy, high aspect ratio and large area structures. In most cases they are fabricated with electron beam lithography (EBL) based techniques plus plasma or wet etching, including extensive efforts in development of the complete nanofabrication process. For this reason, and because of the fact that many devices are made on membranes, a direct patterning technique like FIB offers significant benefits. Here we report on a FZP fabricated in a gold film (Figure 7). It is very difficult to achieve small feature size or high aspect ratio structures on such material, due to gold’s high sputter yield and extreme sensitivity to collateral

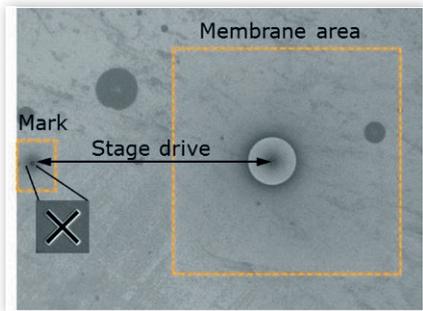


Figure 7: Overview SEM image of the silicon nitride membrane with active area including gold zone plate and reference mark on bulk sample for automatic positioning correction [3]

beam damage. The pattern was fabricated by applying automated position correction with a reference mark on the bulk material far outside the field of view and the sensitive membrane area (Figure 7). This procedure significantly minimized the risk of membrane rupture. The achieved total accuracy of better than 20 nm over 15 h thus includes several highly precise and “blind” drives of the laser interferometer stage to the alignment mark and back to the membrane.

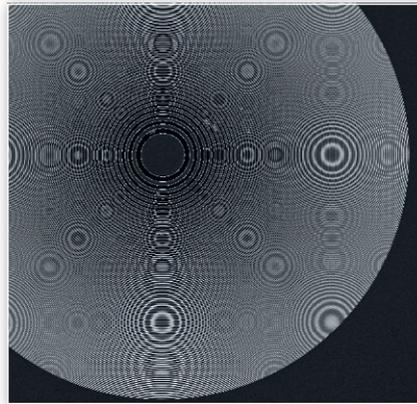


Figure 8: SEM viewgraph of the FZP [3]

Given the dimensions of the FZP, two parameters are the most crucial to investigate: the aspect ratio of the smallest gold zones, and the circularity of the complete zone plate. The first one was checked by cross-sectioning and inspection at four different locations (top, bottom, left, right). The literature shows that an aspect ratio of even 5:1 is an excellent result for patterning of gold. Here the height of the outermost 100 nm zone was found to be almost unchanged (Figure 9).

Verifying the circularity of 100 μm to accuracy of much better than 100 nm is only possible by using a laser interferometer stage in combination with a navigation and metrology software.

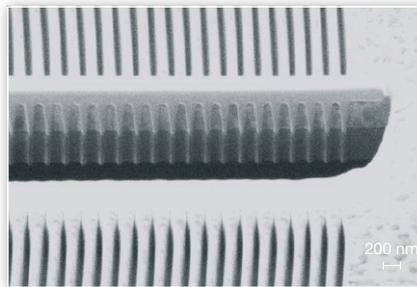


Figure 9: Ion beam cross-section image of outer zones after deposition of a protective and contrast enhancing layer of tungsten [3]

Making use of SEM based metrology, we set up high resolution e-beam line scans at 8 locations including diagonal positions (see Figure 10). We confirmed that the circularity differs by less than 20 nm from ideal, sufficient for excellent FZP performance. This has been confirmed by X-ray tests showing even better resolution than predicted from theory.

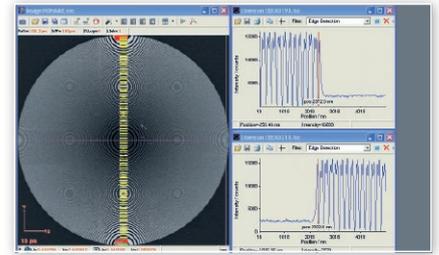


Figure 10: Example of e-beam line scan measurements to determine the diameter in a vertical direction [3]

Combined EBL – FIB process and quality assurance at wafer scale

Membrane based solid state nanopore devices for DNA sequencing and biomolecule analysis are of high interest from both scientific and commercial viewpoints. On the one hand, a feature size below 20 nm for the pore itself is mandatory; on the other, a large number of devices are used for research and, in particular, for commercial screening applications, which requires small batch true wafer scale production.

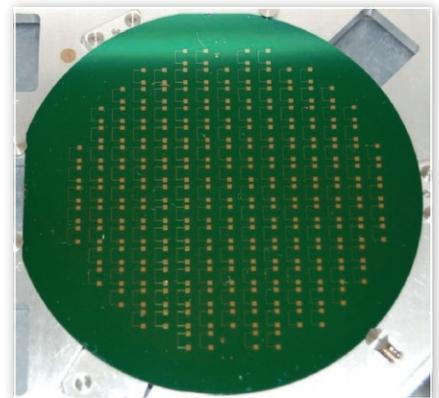


Figure 11: 4 inch wafer with 175 devices of connection pads and read-out electrodes [4]

Here we used a combination of EBL for creating dedicated read-out electrodes for 175 devices on a 4 inch wafer (Figure 11), and FIB for drilling the nanopore well aligned within the corresponding gap (Figure 12). The combination of both techniques, particularly for wafer scale step and repeat and highest resolution processing, is the only solution, since keeping each and every device perfectly in focus can be ensured by the automated height sensing function

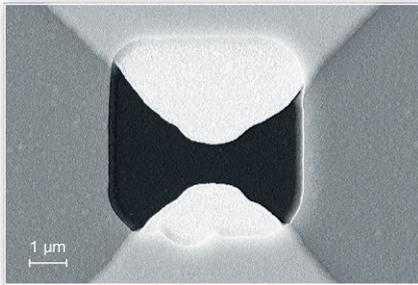


Figure 12: Read-out electrodes within window for placing the nanopore [4]

in Raith systems. Therefore, we were able to achieve 20 nm pores in a 100 nm thick Si_3N_4 membrane across the wafer in a fully unattended way (Figure 13). This has been documented by electron beam quality assurance with a Raith EBL system, which again is possible at wafer scale by means of automation.

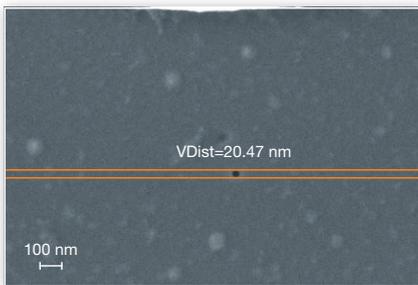


Figure 13: E-beam quality assurance of 20 nm nanopore in 100 nm thick Si_3N_4 membrane [4]

True 3D nanostructures at once over mm length

Finally, structures for micro- and nanofluidic devices are good examples of cases where lateral and depth definition is required to high precision levels and over extended areas. The process for creating a 3D microfluidic mixer channel of constant cross-sectional area while alternating between narrow and deep as well as shallow and wide (Figure 14) can be mastered in a unique approach using FIB nanofabrication. The design elements define the lateral shape and simultaneously the dose for the direct milling process controls the intentionally changing channel depth.

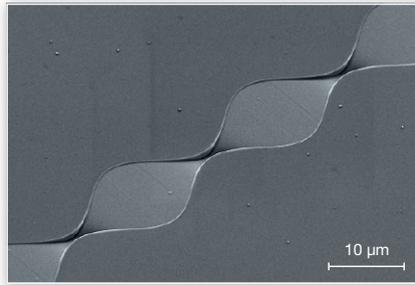


Figure 14: Part of a 1 mm long micro-fluidic mixer channel created by direct milling and stitching [5]

Moreover, employing the laser interferometer stage for stitching multiple base units enables a 1-mm-long channel to be achieved for real devices. This FIB milling process requires some extra care, given the enhanced sputter rate at the edge of subsequent write fields. Possible solutions include minimizing this characteristic by overlapping parts of the write fields with adapted dose distribution (Figure 15). Thus, a very smooth and 3D well defined channel is easily obtained directly in the substrate material.

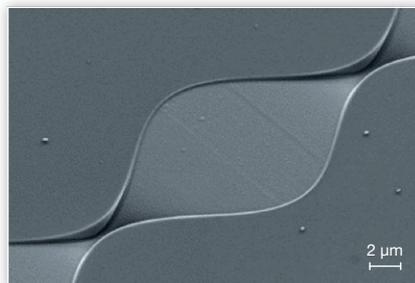


Figure 15: Advanced dose control and overlapping base units during stitching help overcome milling artifacts of adjacent write fields [5]

References

Raith is happy to acknowledge its collaboration with the following partners:

- [1] Trinity College, Dublin
- [2] University of Stuttgart
- [3] Max-Planck-Institute, Stuttgart
- [4] Imperial College London
- [5] Argonne National Laboratory, Chicago

Final Note

Raith's FIB-SEM systems are designed as the first FIB-centric nanofabrication tools for uncompromised state-of-the-art nanopatterning performance. A new world of innovative nanofabrication is now accessible:

- Simplifying process workflow requiring fewer steps, using a large variety of direct FIB patterning techniques
- Highest resolution Ga FIB milling and densely packed nanostructures possible by a low beam tail FIB profile
- Fast FIB fabrication employing ion beam implantation and functionalization
- Creating direct 3D structures in functional layers and resistless patterning on complex or highly topographic devices
- Multiple ion species technology for non-Ga FIB capabilities
- Realizing complex GDSII designs and large area patterns with FIB, including write field stitching and unique continuous patterning modes
- Long-term processes with advanced system stability, enabling the use of smaller beam currents for further improved resolutions
- Lithography class automation including focus control for small batch nanofabrication and unattended FIB jobs on wafer scale
- *In-situ* process control by FIB cross sectioning and SEM inspection
- State-of-the-art EBL patterning and integrated combination of EBL and FIB steps

Explore the enormous potential of Raith's FIB-SEM instruments. Get back to Raith to learn more about the product and its application benefits and to discuss appropriate solutions for your research and development tasks.



Sales

Head office

Raith GmbH
Konrad-Adenauer-Allee 8
44263 Dortmund, Germany
Phone +49 231 95004 0
Email sales@raith.com

Support Europe/Rest of world

Phone +49 231 95004 499
Email support@raith.com

America

Raith America Inc., Islandia, NY
Phone +1 631 738 9500
Email sales@raithamerica.com

Support America

Phone +1 631 738 9500
Email support@raithamerica.com

Asia / Pacific

Raith Asia Ltd., Hong Kong
Phone +852 2887 6828
Email sales@raithasia.com

Support Asia / Pacific

Phone +852 2887 6828
Email support@raithasia.com

China

Raith China Co., Ltd., Beijing
Phone +86 10 828679 22
Email sales@raithchina.com

Support China

Phone +86 10 828679 22
Email support@raithchina.com

India

Raith India Pvt. Ltd., Bangalore
Phone +91 80 2838 4949
Email sales@raithindia.com

Support India

Phone +91 80 2838 4949
Email support@raithindia.com

For further contact information,
please visit our website:

www.raith.com

02-2018

Your challenge is our mission.

RAITH
NANOFABRICATION